


MULTICOM

Clarity Hardware ENG

Code/Rev.: M110/100A Date: 2025-10-22

Phone: +420 251 013 400 clarity@dataapex.com www.dataapex.com DataApex Ltd.
Petrzilkova 2583/13
158 00 Prague 5
Czech Republic

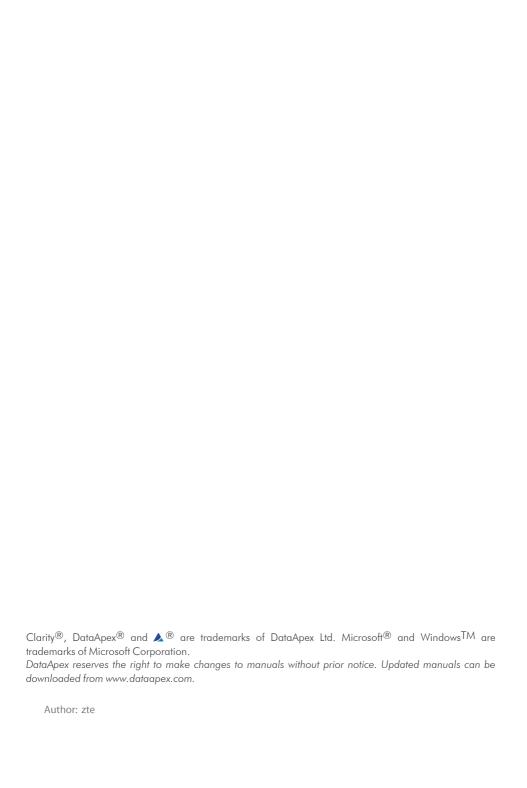

MultiCOM Table of Contents

Contents

1 MultiCOM converter	1
1.1 Side panel	2
1.2 Brief technical overview	2
2 Requirements	3
3 Installation	4
3.1 The MultiCOM Converter with Clarity	4
3.2 The MultiCOM Converter without installing Clarity	5
4 Troubleshooting	6
4.1 Reinstalling - MultiCOM	6
4.2 Changing COM port assignment	10
4.3 How to clean up COM ports that are "(in use)"	12
5 Tables and specifications	15
5.1 Description of the DB9M connector	15
5.2 Disposal and Recycling	16

MultiCOM Table of Contents

To facilitate the orientation in the **MultiCOM** manual and **Clarity** chromatography station, different fonts are used throughout the manual. Meanings of these fonts are:

Open File (italics) describes the commands and names of fields in **Clarity**, parameters that can be entered into them or a window or dialog name.

WORK1 (capitals) indicates the name of the file and/or directory.

ACTIVE (capital italics) marks the state of the station or its part.

Chromatogram (blue underlined) marks clickable links referring to related chapters.

The bold text is sometimes also used for important parts of the text and the name of the **Clarity** station. Moreover, some sections are written in format other than normal text. These sections are formatted as follows:

Note: Notifies the reader of relevant information.

Caution: Warns the user of possibly dangerous or very important information.

Marks the problem statement or trouble question.

Description: Presents more detailed information on the problem, describes its causes,

etc

Solution: Marks the response to the question, presents a procedure how to remove it.

MultiCOM 1 MultiCOM converter

1 MultiCOM converter

This manual describes the use of the MultiCOM USB to RS232 converter.

MultiCOM is a device which connects to the PC via the USB port and provides 6 serial 9-pin ports to the computer. It is also equiped with a free USB port for the connection of the HW key.

MultiCOM converter has been developed for the handling of devices in the LC and GC systems controlled via the RS232 serial interface. When connected to a standard PC, the **MultiCOM** needs no power supply (it is powered from the **USB**), but can also be powered (if needed) from the external power supply adapter.



Fig. 1: MultiCOM - face

The side panel contains two **USB** connectors (one for the connection to the PC, the other for the **HW** key) and the power supply connector. The front panel contains six standard serial 9-pin (plug) ports, each with its own LED communication state indicator.

MultiCOM 1 MultiCOM converter

1.1 Side panel



Fig. 2: Side panel

Green LED status:

- No light not connected to USB, the driver is not installed or is in suspend mode.
- Continual light idle state, no communication.
- Blinking 10100000 (two short consecutive blinks) only sending data (from the USB to the COM port).
- Blinking 01011111 (two short consecutive drop-outs in the continual shine) –
 only receiving data (from the COM port to the USB).
- Blinking 1:1 (2 Hz frequency) both sides are in communication, receiving and sending data.

1.2 Brief technical overview

- 6 RS-232 (non-isolated fully configurable serial ports DB9M).
- 1 USB port for the HW key (or another device with low power consumption).
- 1 USB port for connection to the PC.
- · Powered from **USB** or external power supply.
- USB ports has USB-A connectors and are fully compliant with USB Spec Rev 1.1.

MultiCOM 2 Requirements

2 Requirements

 The MultiCOM converter can be used on any PC with the following operating systems: MS Windows 11/10/8.1/7/Vista/XP/2000 (incl. the 64-bit Windows 11, 10, 8.1, 7).

• The PC must have a free **USB** port (either USB 1 or USB 2 port).

Caution: It is highly recommended to connect the MultiCOM converter directly to the PC (not via USB hubs etc.)

MultiCOM 3 Installation

3 Installation

Ensure that you have **Administrator** access rights in your Windows OS before you proceed with the installation.

Caution: It is necessary to run the pre-installer before the installation (at the time of the first connection of the **MultiCOM** device) to the computer.

There are two possibilities how to install the **MultiCOM** pre-installer into the PC:

- · as a part of the Clarity installation or
- separately as a standalone product

Note: Created COM ports are numbered from the first unused COM number (usually COM3).

3.1 The MultiCOM Converter with Clarity

· Install Clarity from the Clarity installation media.

Clarity ver. 8.2 and newer:

- MultiCOM driver is part of the Typical and Full installation. In case you are installing Custom installation in the Choose Component dialog, unfold the Utils section and check the Multicom driver.
- Continue with the installation process.

Clarity from ver. 2.7 to 8.1:

- Either choose the FULL installation which installs appropriate drivers by default or on the Choose Components screen unfold the Utils section and check the Multicom driver.
- Continue with the installation process.

Clarity ver. 2.6 and older:

- When selecting the type of installation, choose Custom.
- On the next screen, the Utils section (on the bottom of the list) is checked and grayed. Select it and click the Details... button.
- In the available selection, check the checkbox next to the MultiCOM drivers item and press the OK button. Follow the installation process to its end.
- Connect the MultiCOM device to the PC via the supplied USB cable. "Found New Hardware" message will appear on the Windows taskbar and the LED diodes on the MultiCOM will slowly start to light up. Wait until all shine and the taskbar displays the message "New hardware is installed and ready to use".

MultiCOM 3 Installation

3.2 The MultiCOM Converter without installing Clarity

 Install the pre-installer found on the MultiCOM installation media (run the DPINST.EXE file).

Caution: Install the pre-installer before plugging in the MultiCOM device.

Note: The MultiCOM pre-installer on the Clarity installation media can only be used during the installation of the Clarity software. If new installation/re-installation is needed, the MultiCOM installation media should be used instead or the driver

 Connect the MultiCOM device to the PC via the supplied USB cable. "Found New Hardware" message will appear on the Windows taskbar and the LED diodes on the MultiCOM will slowly start to light up. Wait until all shine and the taskbar displays the message "New hardware is installed and ready to use".

should be downloaded from the **DataApex** web pages www.dataapex.com.

4 Troubleshooting

Unplugging the MultiCOM device from the PC while the PC runs may cause the computer to crash in Windows 2000 operating system with the Service Pack installed.

Solution: Unplug the MultiCOM from the USB only after the PC has been turned off.

4.1 Reinstalling - MultiCOM

Normally the drivers are installed with the **FULL** installation option of **Clarity**. In case they have not been installed, **MultiCOM** will not function properly.

"Windows found driver for your device but encountered an error while attempting to install it. The system cannot find the path specified."

Description: This error message occurs when drivers for MultiCOM device have not been

installed and will not function properly.

Solution: Unplug the **MultiCOM** and plug in again. If this does not solve the problem follow the steps below to reinstall/update the **MultiCOM** drivers:

 Locate the item in *Device Manager*. Notice that drivers which are not correctly installed have a warning sign next to them as seen in Fig. 3 on pg. 6.

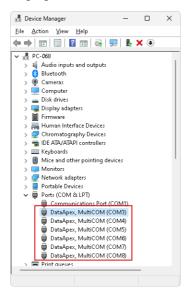


Fig. 3: Device Manager Window

 Right mouse click the item (e.g. DataApex, MultiCOM (COM11)) and select Properties. Dialog showing details about the driver will be opened. On the Driver

tab, select Update Driver...

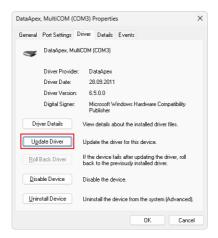


Fig. 4: Dialog for driver properties

• Select "Browse my computer for driver software" option.

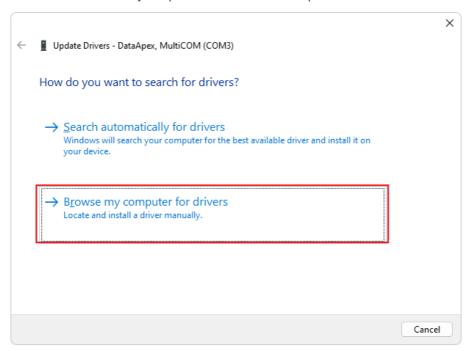


Fig. 5: Dialog for browsing drivers

• Select "Let me pick from a list of device drivers on my computer" option.

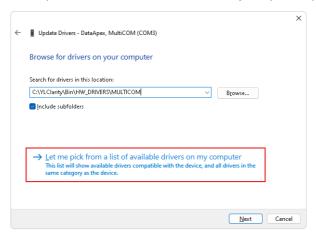


Fig. 6: Dialog for driver properties

· Select "Have Disk" option.

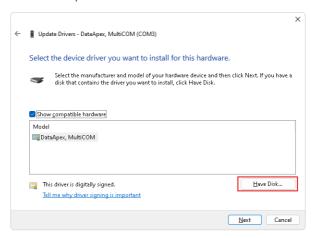


Fig. 7: Dialog for selecting device driver

 Locate the driver in the installation directory, HW_DRIVERS\MULTICOM subfolder and press 2x Open.

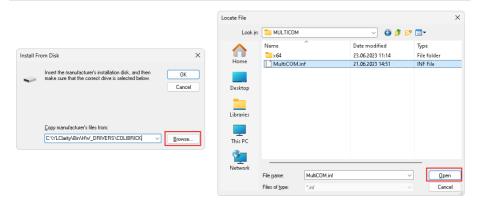


Fig. 8: Dialog for locating the driver Setup Information filetype

• Finish the reinstallation by pressing the *Next* button

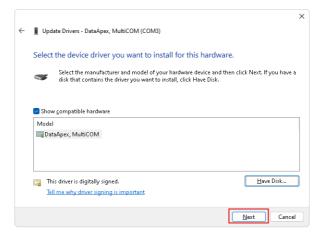


Fig. 9: Dialog for finishing driver reinstallation.

 Upon successful reinstallation of drivers for MultiCOM you will be presented with dialog as seen in Fig. 10 on pg. 10.

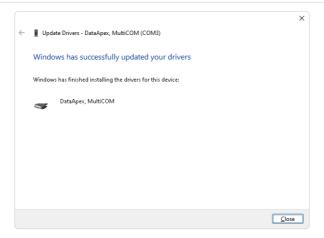


Fig. 10: Successful reinstallation of a MultiCOM driver.

Repeat the aforementioned steps for any other non-functional drivers.

Caution Reinstall/update of drivers may result in assigning a different COM number, for correct communication with Clarity check the appropriate setup.

 If you encounter any difficulties, please contact DataApex Support (www.dataapex.com).

4.2 Changing COM port assignment

In case there is a need to change the **COM** port assignment of the **MultiCOM** device, it can be done via the *Device Manager*.

Changing the assigned COM port number.

Description: In some cases when a new USB device is plugged into the USB port a new COM port is assigned to it. Some drivers do not support COM port numbers higher than 100.

Solution: To change the assigned **COM** port follow the steps bellow:

Go to Device Manager and locate the item where you want to change the
assigned COM port. Right mouse click the item (e.g. DataApex, MultiCOM
(COM8)) and select Properties. Dialog with details about the driver will be
opened. On the Port Settings tab click on Advanced button.

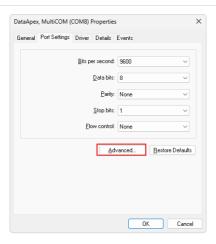


Fig. 11: Device Manager - Port Settings tab

Fig. 12 on pg. 11. displays the Advanced Settings for specific COM port. To change COM port from the default, choose a COM port at ② using the drop-down list.

Note

Already assigned **COM** ports have next to the **COM** port number inscribed "(in use)". Assigning **COM** port that is already "(in use)" may result in malfunction and will trigger a message as seen in **Fig. 13** on pg. **12**.. Clicking **Yes** will change the **COM** port to the selected one from the drop-down list. It is advised to assign only free **COM** ports i.e. **COM** ports not followed by "(in use)".



Fig. 12: Advanced Settings

 Clicking OK will change the assignment of the selected COM port to the selected one from the drop-down list.

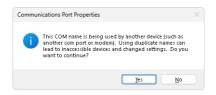


Fig. 13: Message of assigning already used COM port

- Once the COM port has been successfully changed you can locate the new assigned COM port in the Device Manager.
- To change the assignment of other COM ports, repeat previous steps.

4.3 How to clean up COM ports that are "(in use)"

Over time it may happen that excessively large number of COM ports is "in use".

Cleaning up the COM ports that are in "(in use)".

Description: When COM port number is above 100, e.g. COM120, it may happen that the communication between MultiCOM device and Clarity will not work. This is caused by the fact that most control modules do not support COM port higher than

100.

Solution: Before manually assigning the COM ports as described in the chapter "Changing COM port assignment" on pg. 10. You will first need to clean up the COM ports " (in use)". Follow the steps below:

In the Start menu, type regedit and hit enter. You will be prompted with a
message: "Do you want to allow the following program to make changes to this
computer?" After clicking Yes, window as seen in Fig. 14 on pg. 12. will be
opened.

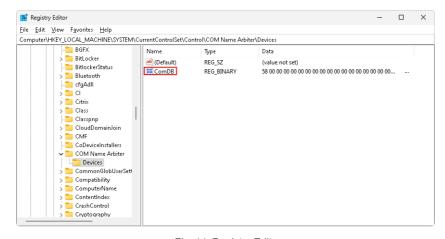


Fig. 14: Registry Editor

- Navigate to Computer\HKEY_LOCAL_MACHINE\SYSTEM\
 CurrentControlSet\Control\COM Name Arbiter
- Right mouse click on the ComDB name and select Modify.... You will be presented with similar data as in Fig. 15 on pg. 13.

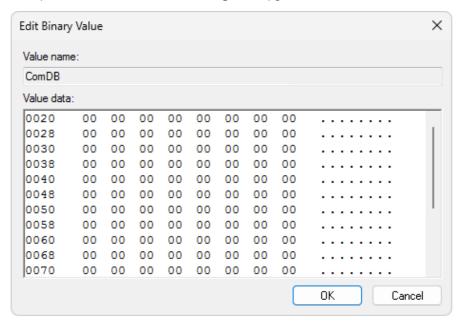


Fig. 15: Registry Editor - modifying ComDB

• Select all the data in the field Value data and delete them manually. Click OK. When you try to Modify... the ComDB again you should be presented with cleaned up window as seen in Fig. 16 on pg. 13..

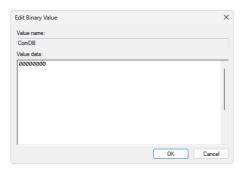


Fig. 16: Registry Editor - cleaned up ComDB

• After cleaning up the **COM** ports you need to restart your **PC**.

- After the restart all the **COM** ports should be free.
- To assign a different COM port, see the chapter "Changing COM port assignment" on pg. 10.

5 Tables and specifications

5.1 Description of the DB9M connector

Description of the DB9M connector

The particular pins of the MultiCOM DB9M connectors are used as follows:

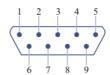


Fig. 17: DB9M connector scheme

Tab. 1: Description of the connector pins:

Pin	Name	Direction	Description
1	CD	←	Carrier Detect
2	RXD	←	Receive Data
3	TXD	\rightarrow	Transmit Data
4	DTR	\rightarrow	Data Terminal Ready
5	GND		System Ground
6	DSR	\rightarrow	Data Set Ready
7	RTS	←	Request to Send
8	CTS	←	Clear to Send
9	RI	←	Ring Indicator

Tab. 2: Complete possibilities of channel configuration:

Parameter	Value		
Data Bits	5, 6, 7 or 8		
Stop Bits	1, 1.5 ^a or 2		
Parity Type	None, Even, Odd, Mark, Space		
Baud Rates	300, 600, 1200, 1800, 2400, 4000, 4800, 7200, 9600, 14400, 16000, 19200, 28800, 38400, 51200, 56000, 57600, 64000, 76800, 115200, 128000, 153600, 230400, 250000, 256000, 460800, 500000, 576000, 921600 ^b		

Notes:

a - 5-bit only

b – 7 or 8 data bits only

Tab. 3: Technical data:

Parameter	Value	
Converter type:	USB to RS232	

Parameter	Value
Number of COM ports:	6, unisolated
USB ports:	1 port to the PC, 1 free port for the HW key or other low consumption device
USB port connectors:	USB-A
Dimensions:	120 x 105 x 22 mm
Weight:	280 g
Power supply:	from the PC via the USB cable or external: stabilized 6V/0.5A
External power supply connector:	plug 1.3mm (outer diameter 3.4 or 3.5 mm)

5.2 Disposal and Recycling



Electrical and electronic equipment must **not be** disposed of as unsorted municipal waste.

It should be collected separately to ensure proper treatment and recycling. Please return used equipment to designated collection points.