


# **TELEDYNE FOXY ISCO 200**

Clarity Control Module ENG

Code/Rev.: M165/100A Date: 2025-10-22

Phone: +420 251 013 400 clarity@dataapex.com www.dataapex.com DataApex Ltd.
Petrzilkova 2583/13
158 00 Prague 5
Czech Republic

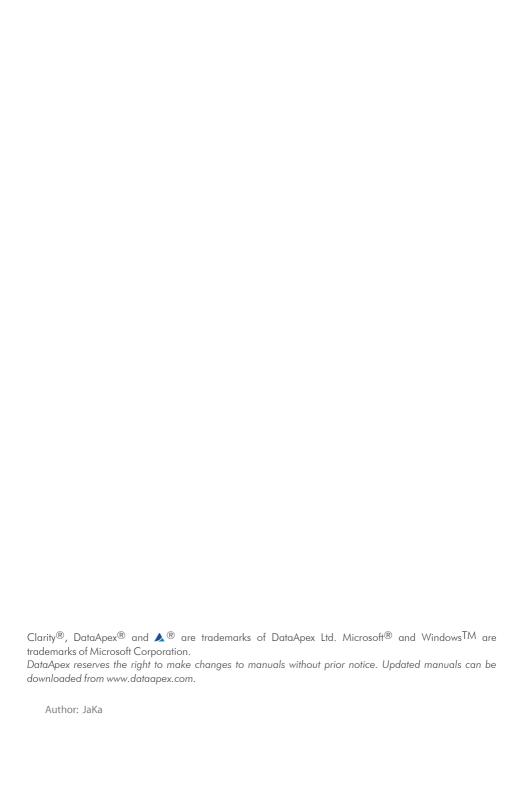

# **Contents**

1 Teledyne Foxy ISCO 200 Control Module	1
2 Requirements	2
3 Installation Procedure	3
3.1 Teledyne ISCO Foxy 200 communication	3
3.2 Clarity Configuration	5
4 Using the control module	8
4.1 Method Setup - FC - Fraction Table	9
4.2 Method Setup - FC - Vial Numbers	13
4.3 Hardware Configuration	15
4.4 Device Monitor	16
4.5 Data Acquisition window	17
4.6 DataApex UNI Setup	18
5 Report Setup	20
6 Troubleshooting	21

To facilitate the orientation in the **Teledyne Foxy ISCO 200** manual and **Clarity** chromatography station, different fonts are used throughout the manual. Meanings of these fonts are:

Open File (italics) describes the commands and names of fields in **Clarity**, parameters that can be entered into them or a window or dialog name.

WORK1 (capitals) indicates the name of the file and/or directory.

ACTIVE (capital italics) marks the state of the station or its part.

Chromatogram (blue underlined) marks clickable links referring to related chapters.

The bold text is sometimes also used for important parts of the text and the name of the **Clarity** station. Moreover, some sections are written in format other than normal text. These sections are formatted as follows:

Note: Notifies the reader of relevant information.

Caution: Warns the user of possibly dangerous or very important information.

#### Marks the problem statement or trouble question.

Description: Presents more detailed information on the problem, describes its causes,

etc

Solution: Marks the response to the question, presents a procedure how to remove it.

# 1 Teledyne Foxy ISCO 200 Control Module

This manual describes the setting of the **Teledyne Foxy ISCO 200** fraction collector. The control module enables direct control of the instrument over serial line.



Fig. 1: Teledyne ISCO Foxy 200 fraction collector

Direct control means that the fraction collector can be completely controlled from the **Clarity** environment. The instrument method controlling the analysis conditions will be saved in the measured chromatograms.

The control is performed via the **UNI Ruby** control module and the **Teledyne Foxy ISCO 200** script.

# 2 Requirements

- Clarity Installation USB with LC control (p/n A24) license allowed.
- Free serial COM port in the PC.

Note: Mode

Modern computers usually have only one (if any) serial (COM) port installed. To use more devices requiring the RS232 port, the **MultiCOM** adapter (p/n MC01) is available.

Serial DB9F-DB25M cable (p/n SK12).

Note: Cables are not part of Clarity Control Module. It is strongly recommended that you order the required cables together with the Control Module.

 Firmware version of the Teledyne Foxy ISCO 200 collector must be either 2135-061 (standard version) or 2135-063 (CombiFlash version).

Note: Other firmware versions may function but have not been tested by **DataApex**.

# **3 Installation Procedure**

# 3.1 Teledyne ISCO Foxy 200 communication

The **Teledyne Foxy ISCO 200** is controlled by serial (RS232) communication. It uses a serial DB9F-DB25M straight cable (p/n SK12) described in the picture below.

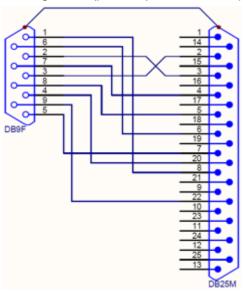


Fig. 2: Serial straight DB9F-DB25M cable

On top of that, the fraction collector must be switched to the same Baud Rate as set in **Clarity** and the ID=5 to work with the control module. To do so, start the fraction collector and load the *Program A*, then select the *Edit A* option. Select page 3 from the Edit submenu and set the *Serial* function, then set the Baud Rate as desired and *ID* to 5 in the next screen.

Caution: The fraction collector must be set to precise settings, refer to procedure given below.

- · Switch ON the fraction collector.
- Press PROGRAM button.
- Press A button in order to edit Program A.
- Press A button again in order to edit item TIME\DROP\VOL.
- Press C button within next screen (use of time windows/peak detection within program) on order to select No option.
- Press A button to set Rack 1 item.
- Keep pressing A button in order to set Tube Types item. The preset Tube
  Type has to correspond to type of tubes present in Rack 1.
- Press D button to set number of tubes in Rack. Number of tubes should be set using numerical keyboard below display and be confirmed by pressing of ENTER button.
- Set 1:00 and Press ENTER in next Fraction Size screen.
- On the next screen is necessary to set Rack 2 if used. Pressing of button B or C or D has to correspond to reality.
- For the following steps of this procedure it is assumed there is used only Rack
  1 on the fraction collector.
- For the following steps of this procedure is assumed there is used only Rack 1
  on the fraction collector, thus button D should be pressed.
- Keep pressing **D** until there is reached *Page* 3 in Menu.
- Press C button in order to edit SERIAL item.
- Press C button in order to select 9600 item.
- Set 5 value as ID using numerical keyboard below display and confirm by pressing ENTER button.

#### Caution:

In case there is changed type of tubes used (in any rack) it is necessary edit settings accordingly in the <u>Foxy 200 Setup</u> and in the used fraction collector's Program as well.

In case there is changed number of tubes used (in any rack) within *Program* of the fraction collector it is necessary consider this value during editing (item *Last Vial Number*) of the <u>Vial Numbers tab in the Method Setup dialog</u>.

# 3.2 Clarity Configuration

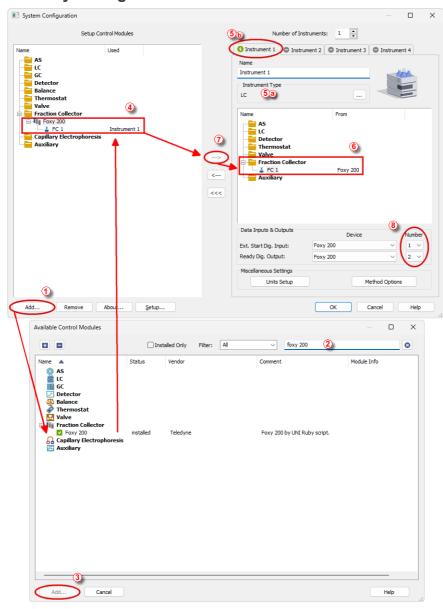


Fig. 3: How to Add Foxy 200 Control Module

- Start the **Clarity** station by clicking on the **\( \rightarrow**\) icon on the desktop.
- Invoke the System Configuration dialog accessible from the Clarity window using the System Configuration... command.
- Press the Add button (1) on 3.2 on pg. 5.) to invoke the Available Control
   Modules dialog.
- You can specify the search filter ② to simplify the finding of the driver.
- Select the correct item and press the Add (③ on Fig. 3 on pg. 5.) button. Each
  device with already created profile should have its own item named accordingly
  in the Available Control Modules dialog.
- The DataApex UNI Setup dialog will appear.

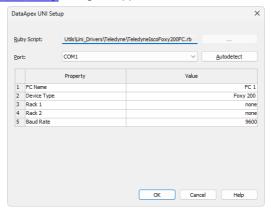


Fig. 4: DataApex UNI Setup - Teledyne ISCO Foxy 200

- Select the appropriate script using the \_\_\_\_\_\_ button. The scripts already prepared in Clarity are stored in the UTILS/UNI\_DRIVERS/ sub-directory (sorted in further sub-directories by the instrument manufacturers) of the Clarity installation directory.
- Set the correct communication *Port* and click on the *AutoDetect* button to establish communication with the device.
- You may fill in the custom *Device Name*.

Note: The DataApex UNI Setup dialog is described in detail in the chapter "DataApex UNI Setup" on pg. 18.

- The **Teledyne Foxy ISCO 200** item ④ will appear in the *Setup Control Modules* list of the *System Configuration* dialog.
- Change the *Instrument Type* 3 on the desired *Instrument* tab 3 to LC.
- Drag the appropriate item from the Setup Control Modules list on the left 4 to the Instrument tab on the right 6 , or use the button 7.

• Set the Ext. Start Dig. Input and Ready Dig. Output numbers <sup>®</sup> for your acquisition card according to the wires used for synchronization.

# 4 Using the control module

After adding and setting up the fraction collector a new tab *Method Setup - FC* will appear in the *Method Setup* dialog. A new **Teledyne Foxy ISCO 200** FC section enabling the monitoring of the current fraction collector state will be also created in the **Device Monitor** window.

## 4.1 Method Setup - FC - Fraction Table

The *Method Setup - FC - Fraction Table* tab is used for setting the automatic fraction collection program of the **Teledyne Foxy ISCO 200** fraction collector. To display the actions of the fraction collector in *Data Acquisition* window or *Chromatogram* window, click the right mouse button on the graph, choose *Properties...* to open the **Graph properties** dialog and check the *Show Events* option.

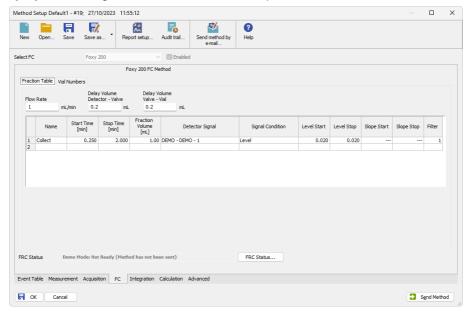


Fig. 5: Method Setup - FC - Fraction Table

#### Flow Rate

Specifies the flow rate of the liquid going to the fraction collector. Together with *Fraction Volume* parameter from the *Fraction Table* it is used to prevent vial overflow.

### **Delay Volume Detector - Valve**

Defines the volume of tubing between the detector and the fraction collector's waste/collect valve.

The delayed volume is displayed in the <u>Data Acquisition</u> window as a vertical line trailing the signal front. Choose *Properties*... to open the *Graph Properties* dialog and check the *Show Events* checkbox.

#### **Delay Volume Valve - Vial**

Defines the tubing volume between the waste/collect valve and the vial in the fraction collector tray. If the fraction collector has no waste/collect valve, leave this field empty and enter the total tubing volume from the detector and fraction collector tray in the Delay Volume Detector - Valve field.

#### Fraction Table

Sets up the fraction collector behavior. The table rows run in chronological order and time intervals should not overlap. If a collection is active when the interval ends, it stops immediately. Collection based on negative peaks is not supported; to collect such peaks, invert the detector signal before setting up the method.

#### Name

This name will be displayed in the *Special Results* table for all fractions collected in the given interval.

#### Start Time [min]

Beginning of the time interval where the Signal Condition will be evaluated.

#### Stop Time [min]

End of the time interval where the *Signal Condition* will be evaluated.

#### **Fraction Volume**

Defines the volume assigned to a single fraction. The value should not exceed the physical vial capacity to avoid overflow. When the collected volume reaches the set value, the fraction collector automatically switches to the next vial.

#### **Detector Signal**

Specifies the detector signal used to trigger fraction collection according to the *Signal Condition* in the *Fraction Table*. Each row of the table can be linked to a different detector signal. Only detector signals that are enabled for acquisition are available for use in the *Fraction Table*.

#### **Signal Condition**

Signal Condition is evaluated within the time interval defined for each row in the Fraction Table. If the condition is met multiple times during the interval, fraction collection starts each time, with the vial automatically changed after every stop. If the condition is not met at all, no collection is performed for that interval.

None - The fractions are collected for the entire time interval defined by Start Time and Stop Time regardless of the detector signal.

Note:

If two Fraction Table rows follow each other directly in time (Stop Time = Start Time) and both have the Signal Condition = None, the FC does not switch to Waste between the rows. Skipping the switch to Waste is possible only if the collector still has a free vial available and the Waste During option in the Vial Change method is not selected.

#### Level - The fraction collection:

- starts when the detector signal increases above the Level Start
- stops when the detector signal decreases below the Level Stop

#### Slope - The fraction collection:

- starts when the first derivative of the detector signal increases above the Slope Start
- stops when the first derivative of the detector signal first decreases and then increases above the negative value of the Slope Stop. The following image

displays the course of the first derivative and demonstrates *Slope Stop* activation. Carefully read the slope values at all *End Fraction* points and observe the course of the first derivative before it reaches these points.

For accuracy, the *Slope Stop* condition is applied with a tolerance of 5%. The entered value is multiplied by 1.05 (e.g., an entered value of 0.100 is evaluated as 0.105).

It is strongly recommended **not** to use *Slope Signal Condition* for fraction collection based on noisy detector signal.

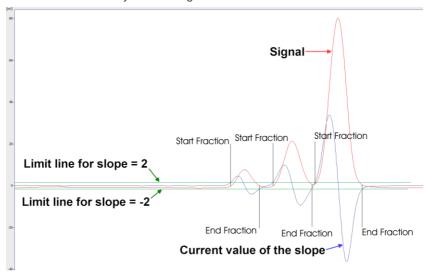


Fig. 6: First derivative course

#### Start L+S End L+S - The fraction collection:

- starts when both start signal conditions are fulfilled simultaneously, detector signal increases above Level Start value and also the first derivative of the detector signal increases above Slope Start value.
- stops when both stop signal conditions are fulfilled simultaneously, detector signal decreases below Level Stop value and also the first derivative of the detector signal first decreases and then increases above the negative of the Slope Stop value. If only one of the stop signal conditions is fulfilled the fraction collection will not stop until Stop Time is reached.

#### Start L+S End L or S - The fraction collection:

 starts when both start signal conditions are fulfilled simultaneously, detector signal increases above Level Start value and also the first derivative of the detector signal increases above Slope Start value.  stops when any of the stop signal conditions is fulfilled (whatever occurs first), detector signal decreases below *Level Stop* value or the first derivative of the detector signal first decreases and then increases above the negative of the *Slope Stop* value. If none of the stop signal conditions is fulfilled the fraction collection will not stop until *Stop Time* is reached.

#### Start L or S End L+S - The fraction collection:

- starts when any of the start signal conditions is fulfilled (whatever occurs first), detector signal increases above Level Start value or the first derivative of the detector signal increases above Slope Start value.
- stops when both stop signal conditions are fulfilled simultaneously, detector signal decreases below Level Stop value and also the first derivative of the detector signal first decreases and then increases above the negative of the Slope Stop value. If only one of the stop signal conditions is fulfilled the fraction collection does not stop until Stop Time is reached.

#### Start L or S End L or S - The fraction collection:

- starts when any of the start signal conditions is fulfilled (whatever occurs first), detector signal increases above Level Start value or the first derivative of the detector signal increases above Slope Start value.
- stops when any of the stop signal conditions is fulfilled (whatever occurs first), detector signal decreases below *Level Stop* value or the first derivative of the detector signal first decreases and then increases above the negative of the *Slope Stop* value. If none of the stop signal conditions is fulfilled the fraction collection does not stop until *Stop Time* is reached.

#### Filter

Sets the data filter for the detector signal defined in the corresponding row of the *Fraction Table*. The default value of 1 means no filtering, value of X means that every X incoming data points are averaged into one point used for evaluating fraction start/stop conditions. The higher the *Filter* value, the greater the delay in responding to signal conditions, especially with detectors that have a low data rate.

Filter settings influence the value that should be set to the *Slope Start* and *Slope Stop* columns. To determine the correct trigger values for the *Slope Start* or *Slope Stop* with the *Show Slope/Level* tool in the *Chromatogram* window, the *Global Filter – Bunching* value in the *Integration Table* should match the *Filter* value set here.

The values coming from *Detector Signal* are evaluated in bunch approximately every 300 milliseconds with respect to Start and Stop values inserted in the *Fraction Table*. As a result, the positions of fraction collection start and stop events may vary by  $\pm 300$  ms between chromatograms with identical signal curves. For detectors with a high sampling rate (e.g., 50 Hz), this corresponds to a variation of up to  $\pm 15$  data points (50 Hz = 50 points per second; 300 ms  $\approx 15$  points).

## 4.2 Method Setup - FC - Vial Numbers

The *Method Setup - FC - Vial Numbers* tab is used for setting the behavior of the **Teledyne Foxy ISCO 200** fraction collector.

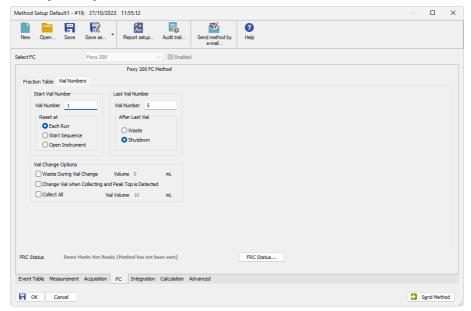


Fig. 7: Method Setup - FC - Vial Numbers

#### Start Vial Number

#### Vial Number

Sets the number of the vial where the first collected fraction will be stored. The current vial number is incremented each time the *Next* signal is activated. The vial number is stored in the *Results Table*.

Note:

The **Teledyne Foxy ISCO 200** fraction collector does not allow to move the collection vessel to a specific vial. The collection will always start on the first vial or last vial +1 from the previous run, depending on the *Reset at* settings.

#### Reset at

Specifies the conditions of the current vial number setting to *Start Vial Number* in three different situations:

- Each Run the vial number is reset at the start of each acquisition.
- Start Sequence the vial number is reset only at the start of a sequence.
- Open Instrument the vial number is reset only once, when the Instrument is opened.

#### **Last Vial Number**

#### Vial Number

Sets the number of the last available vial. The **Teledyne Foxy ISCO 200** fraction collector can set the tray used in the <u>DataApex UNI Setup</u> dialog, so the value range of this field is limited.

#### After Last Vial

Specifies the action executed when current vial number exceeds the Last Vial Number

- Waste the acquisition continues, but the fraction table is ignored and the samples are automatically discarded.
- Shutdown the acquisition is immediately stopped and the analysis aborted.

### **Vial Change Options**

### **Waste During Vial Change**

This option prevents spilling when the vial is changed during collection. It temporary stops the fraction collection before changing the vial and restarts it again when the vial is swapped. The vial is changed in the middle of the wasting interval; the volume wasted during vial change can be specified in the *Volume* field. Volume units are the ones set in the <u>Units Setup</u> dialog. This functionality is not applicable (even if checked) when *Collect All* option is selected.

#### Change Vial when Collecting and Peak Top is Detected

If enabled, the vial is automatically swapped when the peak top is detected during collection. Be aware this option is not applied if the active row in the *Fraction Table* is set to *None* in the *Signal Condition* column.

#### Collect All

When checked, fractions are collected during the whole course of the analysis, regardless of the settings in the *Fraction Table*. The FC module then sends the *Next Vial* command **only**, which makes this mode of operation suitable for a fraction collector without a waste position. *Collect All* mode does not permit *Waste During Vial Change*.

The maximum amount of eluent collected in a single vial is limited by the value in *Vial Volume* field. Once this volume is filled, it begins collecting into the next successive vial.

Note:

With Collect All option the valve is not controlled.

Warning:

Once error message "Vial number exceeded last vial number" is displayed, it means that the FC was already at the last position when command Next Vial was received. Collecting stops at position which is equal to the Last Vial Number - 1 because it's the last possible position to collect (due to Waste). Correct appropriately your method.

# 4.3 Hardware Configuration

The *Hardware Configuration* dialog invoked by using the *FRC Status...* button from the *Method Setup - FC* dialog displays the configuration of the **Teledyne Foxy ISCO 200**, namely information about configured racks and communication baud rate.

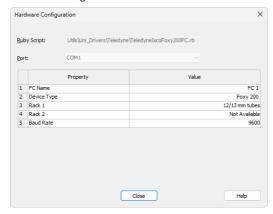


Fig. 8: Hardware Configuration Teledyne ISCO Foxy 200 Fraction Collector

#### 4.4 Device Monitor

The window with the fraction collector status can be invoked by the *Monitor - Device Monitor* command from the *Instrument* window or using the *Operice Monitor* icon.

#### **Device Monitor - Fraction Collector**

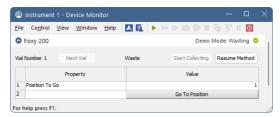


Fig. 9: Device Monitor - Teledyne Foxy 200 Fraction Collector

#### **Next Vial**

Moves the collection vessel one vial forward. Using this button while the collector is performing its program set on the <u>Method Setup - FC - Fraction Table</u> tab during an analysis run immediately overrides the automatic collection program. To return to the automatic program, use the *Resume Method* button.

#### Start Collecting / Stop Collecting

The description on the button changes according to the state in which the fraction collector is - while it is collecting, the *Stop Collecting* button forces the fraction collector to stop collecting, while it is in a waste cycle, the *Start Collecting* button will force the collector to start collecting. Using any of the two actions while the collector is performing the program set on the <a href="Method Setup-FC-Fraction Table">Method Setup-FC-Fraction Table</a> tab during an analysis run, immediately overrides the automatic collection program. To return to the automatic program, use the *Resume Method* button.

#### **Position To Go**

Allows to set the position to which the fraction collector should be switched. The switching itself is performed using the *Go To Position* button.

Note: The button only functions when the fraction collector is not wasting.

#### **Resume Method**

Resumes the actions of the automatic fraction table set on the <u>Method Setup - FC - Fraction Table</u> tab. If the fraction collector is collecting at that moment, it will start a waste cycle and then resume the function of the fraction table.

# 4.5 Data Acquisition window

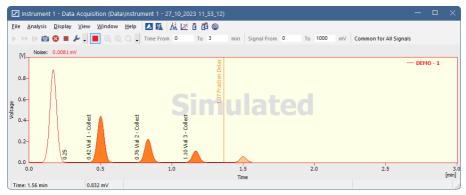


Fig. 10: Data Acquisition window

The *Data Acquisition* window can display the fractions in the graph using the background color and also the start and stop fraction events and marker of the Fraction Delay. To see the fractions and markers right click on the graph, select *Properties*... to open the *Graph Properties* dialog and check the *Show Events* checkbox.

When the Fraction Delay is set, there will be two colors present in the graph - lighter and darker red. The lighter red color shows the expected duration of the fraction (which **Clarity** only presumes based on data points it got so far), while the darker red color shows the placement of the real fractions.

Note: The Fraction Delay in the picture will appear if the Delay Volume field in the Method Setup - FC - Fraction Table dialog is set to a nonzero value.

## 4.6 DataApex UNI Setup

The appearance of the *DataApex UNI Setup* dialog depends on the presence of the selected Ruby Script - if the script is not present, only the *Ruby Script* field is visible.

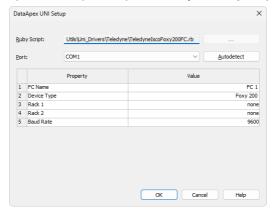


Fig. 11: DataApex UNI Setup - Teledyne ISCO Foxy 200

#### **Ruby Script**

Displays the selected Ruby Script. The correct TELEDYNEISCOFOXY200FC.RB script for the controlled instrument depends on the instrument manufacturer. For Teledyne ISCO instruments there is a subdirectory with the scripts (accessible through the \_\_\_\_\_\_) button of the **Clarity** installation folder (C:\CLARITY\BIN by default). The selection of different script is disabled in the dialog for devices added with pre-selected scripts, to be able to select the script manually, use the **UNI RUBY** item in the auxiliary category.

#### Port

Defines the communication port used, possible values dependent on the type of communication of the device and/or available ports in the PC.

#### AutoDetect

It is used for verifying the device communication over the serial port selected above.

#### **FC Name**

Allows you to set the custom name of the fraction collector. This name (entered into the *Value* column) will be used throughout the **Clarity** station.

#### **Rack Selection**

Allows the user to select the correct rack type installed on the **Teledyne Foxy ISCO 200** fraction collector. **Teledyne Foxy ISCO 200** fraction collector supports two racks. These settings **cannot** be read rom the fraction collector itself by using the *Autodetect* button.

#### **Baud Rate**

Sets the Baud Rate communication parameter for the **Teledyne Foxy ISCO 200** fraction collector. Value set in this dialog has to correspond in method in fraction

### collector itself.

Note:

The setting of the racks to be used is performed in the Method A of the fraction collector and cannot be changed from the side of **Clarity** - this is the limitation of the fraction collector communication protocol.

# **5 Report Setup**

The fraction collector section on the method report can be enabled by checking the *Instrument Control* checkbox on the *Method* tab of the *Report Setup* dialog.

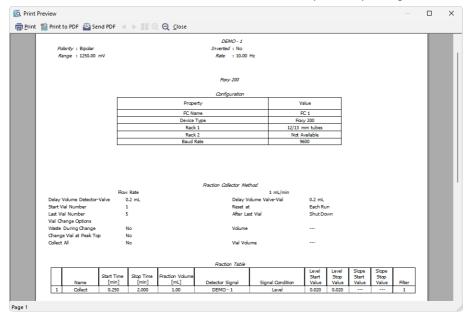


Fig. 12: Report Setup - Teledyne ISCO Foxy 200 Fraction Collector

All of the parameters set in the Method Setup - FC - Fraction Table and Method Setup - FC - Vial Numbers tabs are reported, as well as the custom FC Name and other parameters set in the DataApex UNI Setup dialog.

# 6 Troubleshooting

When the solution to a problem cannot be found easily, a recording of the communication between Clarity and the fraction collector will significantly help DataApex support.

The data recording can be enabled by adding or amending the COMMDRV.INI file in the Clarity installation directory (C:\CLARITY\CFG by default). The file can be edited in any text editor (e.g. Notepad). The following section should be edited or added:

[COM1] echo=on textmode=on filename=Teledyne\_Foxy\_COM1\_%D.txt reset=off

Note: Instead of COM1, type the communication port used to communicate with the fraction collector. This port number is displayed when the FRC Status... button in the Method Setup - FC tab is clicked. The DataApex Uni Setup dialog with correct COM port number is invoked.

> %D (or %d) in the filename parameter means that the log will be created separately for each day. The reset=off parameter disables deleting the content of the log each time the station is started during the same day.

The created \*.TXT files will be of great help in the diagnosis of not documented errors and communication issues.

#### Note: