

U-PAD A/D CONVERTER

Clarity Hardware ENG

Code/Rev.: M059/80F Date: 2025-10-22

Phone: +420 251 013 400 clarity@dataapex.com www.dataapex.com DataApex Ltd.
Petrzilkova 2583/13
158 00 Prague 5
Czech Republic

Sections of the manual connected only to the Clarity Full version are marked with the Vill version.	
Clarity [®] , DataApex [®] and \triangle [®] are trademarks of DataApex Ltd. Microsoft [®] and Windows TM of trademarks of Microsoft Corporation. DataApex reserves the right to make changes to manuals without prior notice. Updated manuals can adownloaded from www.dataapex.com.	
Author: zte	

Contents

1 U-PAD acquisition device	1
1.1 Side panel of the U-PAD device	
1.2 A/D converter types in general	1
2 Requirements	3
3 Installation	4
3.1 The U-PAD	4
3.2 Standard cable for Clarity station	6
3.3 Connection with chromatograph	7
3.3.1 Connection of signal cables	7
3.3.2 Connection of starting cables	8
3.4 Clarity Configuration	9
3.4.1 Measuring on multiple instruments	11
3.4.2 Using multiple U-PAD converters	12
4 Using the U-PAD	13
4.1 DataApex U-PAD1 Setup	13
4.2 Method Setup	16
4.2.1 Method Setup - Acquisition	16
4.2.2 Method Setup - Measurement	18
4.3 Digital Inputs and Outputs	19
4.4 Device Monitor	20
5 Troubleshooting	21
5.1 Locate your problem	21
5.2 Problems with U-PAD	22
5.2.1 How to check the U-PAD driver	22
5.3 Manual installation	24
5.3.1 Installation and reinstallation in Windows XP/Vista/7/8.1	24
5.3.1.1 Reinstallation of drivers using the System Restore Point	26
5.3.2 Reinstallation of drivers in Windows 2000	27
5.3.2.1 Deleting incorrect *.INF files in the system	28
5.3.2.2 Installing the correct driver manually	29
5.4 Data Acquisition - non-functional	32
5.5 Data Acquisition - Simulated	34
6 Tables and specifications	35
6.1 Description of the INT7 connector (Male)	35
6.2 Parameters of digital inputs and outputs	
6.3 U-PAD - CE Conformity Declaration	37

U-PAD A/D Converter Table of Contents

6.4 Technical data	. 3
6.5 Disposal and Recycling	.38

U-PAD A/D Converter Table of Contents

To facilitate the orientation in the **U-PAD A/D Converter** manual and **Clarity** chromatography station, different fonts are used throughout the manual. Meanings of these fonts are:

Open File (italics) describes the commands and names of fields in **Clarity**, parameters that can be entered into them or a window or dialog name.

WORK1 (capitals) indicates the name of the file and/or directory.

ACTIVE (capital italics) marks the state of the station or its part.

Chromatogram (blue underlined) marks clickable links referring to related chapters.

The bold text is sometimes also used for important parts of the text and the name of the **Clarity** station. Moreover, some sections are written in format other than normal text. These sections are formatted as follows:

Note: Notifies the reader of relevant information.

Caution: Warns the user of possibly dangerous or very important information.

Marks the problem statement or trouble question.

Description: Presents more detailed information on the problem, describes its causes,

etc

Solution: Marks the response to the question, presents a procedure how to remove it.

1 U-PAD acquisition device

This manual describes the use of the **U-PAD A/D Converter** with the **Clarity** software ver. **2.7 and later**.

U-PAD is an external measuring unit designed for acquisition of data from any chromatograph for the **Clarity** station. **U-PAD** uses the **USB** communication channel and thus is directly powered from the **PC**. The unit can be connected or disconnected anytime without the necessity of switching the computer off.



Fig. 1: U-PAD

Brief technical overview

- · 2 data acquisition channels jointly isolated from USB ground
- 2 digital inputs
- · 2 digital outputs designed also as relay contacts
- USB interface
- · excellent reliability and temperature stability

1.1 Side panel of the U-PAD device

The side panel contains **USB** connector and **green LED** indicating the status of the device



Fig. 2: Side panel of the U-PAD

Green LED status

- Blinking device is Idle
- Steady light at least one channel is acquiring data

1.2 A/D converter types in general

Prior to collecting and processing the chromatographic signal by a computer, the analog signal of the detector must be first converted to a digital form by an analog/digital (A/D) converter.

In principle there are three types of A/D converters – sampling, integrating and $\Delta\Sigma$ (**Delta-Sigma** - with continuous integration).

The **U-PAD** converter uses Delta-Sigma integration.

Advantages of the U-PAD A/D converter

- No loss of input signal during integration. There are no time delays during which the converter fails to integrate the input signal.
- No continuous servicing of the input analog switches, which would otherwise cause errors (offset, noise).
- Substantially smaller amounts of data entering the computer due to individual data items already representing partial integrals.

U-PAD A/D Converter 2 Requirements

2 Requirements

The **U-PAD** external **USB** A/D **Converter** can be used with the **Clarity** software on a PC with a free **USB** slot and a **32-bit Windows OS**.

Caution: U-PAD A/D Converter is supported up to Windows 10 OS.

Tab. 1: Hardware and Software Compatibility:

	Windows			
Software	11	10 (32-bit only)	8.1 (32-bit only)	7 ^{SP1} (32-bit only)
Clarity	No	Yes (from 6.2)	Yes (from 4.0.3)	Yes
Clarity Lite	No	Yes (from 6.2)	Yes (from 4.0.3)	Yes
CSW32	No	No	No	Yes

Tab. 2: Hardware and Software Compatibility with obsolete operating systems:

	Windows			
Software	XP ^{SP3} (32-bit only)	Vista ^{SP2} (32-bit only)	2000	
Clarity	Yes (until 7.4.1)	Yes (until 7.2)	Yes (until 3.0)	
Clarity Lite	Yes (until 7.4.1)	Yes (until 7.2)	Yes (until 3.0)	
CSW32	Yes	No	Yes	

3 Installation

Ensure that you have *Administrator* access rights in your **Windows OS** before you proceed with the installation.

U-PAD A/D Converter is not part of the **Clarity** *Typical* installation. To install it, select the *Full* installation or the *U-PAD* item in the *Choose Components* dialog during the installation of **Clarity**.

3.1 The U-PAD

· Install the Clarity software from the DVD-ROM.

Caution: Install Clarity before plugging in any devices.

The drivers were copied to your computer during the installation of the Clarity software.

 Plug the U-PAD device into the USB slot and the Found New Hardware Wizard will appear.

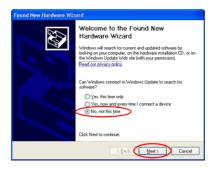


Fig. 3: Found New Hardware Wizard - Step 1

Select No. not this time and click Next.

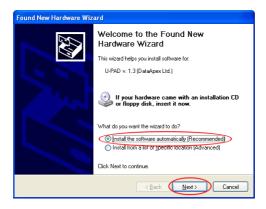


Fig. 4: und New Hardware Wizard - Step 2

• Select Install the software automatically (Recommended).

Note: If you had previously used an older version of the **U-PAD driver**, select Install from a list or specific location (Advanced), in the next step select Don't search, I will choose the driver to install and proceed analogically to re-installation procedure described in the chapter "Installation and reinstallation in Windows

XP/Vista/7/8.1" on pg. 24.

• During the installation process, the *Hardware Installation* dialog with information that the software for **U-PAD** has not passed the Windows Logo testing may appear.



Fig. 5: Windows Logo Testing Message

• Click the Continue Anyway button.

Finally the Found New Hardware Wizard will display the Completing the Found New Hardware Wizard dialog.

- Click the Finish button.
- The driver is now installed, proceed to the chapter "Standard cable for Clarity station" on pq. 6.

Note: When installing multiple **U-PAD**s it is recommended to restart the PC after the installation to let the USB devices set in the order in which they always appear in the lists

3.2 Standard cable for Clarity station



Fig. 6: Cables for the Clarity station package

The standard **Clarity** station package includes a cable composed of signal and starting cables for connecting the **Clarity** station to the chromatograph.

Signal cables

Labeled **DET1** and **DET2**, the cables are supplied as standard without connectors [bare stripped and tinned endings – red (+), white (-) and shielding (analogue ground)].

Starting (marker) cables

Labeled **IN1** and **IN2**, ended with CINCH connector (female). The following cables can be connected to this CINCH connector:

- A cable terminated with <u>free leads</u> [red (+), shielding (digital ground)] for direct connection to the chromatograph.
- A cable terminated with a <u>button</u> for cases where a starting contact is not available and it is necessary to perform a manual start.

Both types of cables are supplied for each starting cable.

Digital output cables

Relay contacts labeled **OUT1R** and **OUT2R**, terminated with free leads. These are used for synchronizing autosamplers without an **AS Control** module in the *ACTIVE* Sequence or for controlling other devices.

At the converter end the cable is always terminated with a DB-37F connector (37-pin female connector).

3.3 Connection with chromatograph

Connect the cables according to one of the following diagrams in the **Fig. 7** on pg. **7**. Use the <u>symmetrical connection</u> only when you are sure that the chromatograph/detector is equipped with a symmetrical output – it is necessary to read through the instructions for the corresponding chromatograph.

All current **DataApex** A/D Converters **INT7**, **INT9**, **U-PAD**, **U-PAD2** and **Net-PAD** use the same standard **INT7** Connector.

Note: A description of the INT7 connector can be found in the chapter "Tables and specifications" on pg. 35.

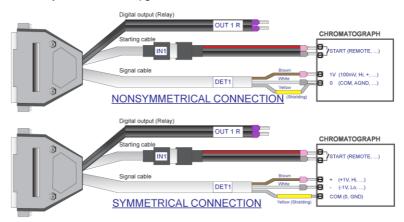


Fig. 7: Connecting Clarity to chromatograph

3.3.1 Connection of signal cables

Signal inputs of the A/D converters are symmetrical: + (red), - (white) and analogue ground (copper braiding).

Caution:

Shielding must be connected. It serves not only as the shielding but also as the analogue earth against which the measurement takes place. In the case of asymmetrical output of a detector (only two connectors) shielding must be connected to the white lead! No lead of the signal cable may remain unconnected.

Try to connect to the detector output of the chromatograph with the largest possible signal level, usually indicated in the chromatograph as **INTEGRATOR** (signal approx. 1 V). The level of the signal on the output marked **RECORDER** is usually only approx. 10 mV. For easier alterations of the wiring, we supply a **SV9 Terminal board** with screw contacts for **INT7**, **INT9**, **U-PAD**, **U-PAD2** and **Net-PAD** A/D converters.

Note:

This **SV9 Terminal Board** is not suitable for applications with small signals or with high electromagnetic interference. This is because the **SV9 Terminal Board** uses leads and screw contacts that are not shielded.

3.3.2 Connection of starting cables

Starting input reacts to a change of the TTL level (5 V / 0 V) or to a connection by any contact (button, contact of relay).

The input implicitly reacts to a change from *HIGH* to *LOW* (contact closure). The input function may be altered by changing the *Down* option to *Up* in the *External Start/Stop* section of the *Method Setup - Measurement* dialog (accessible from the *Instrument* window using the *Method - Measurement* command).

For schemes and hints describing the typical autosampler wiring, please see the **Connecting Autosamplers** chapter of the **Getting Started Manual**.

3.4 Clarity Configuration ______

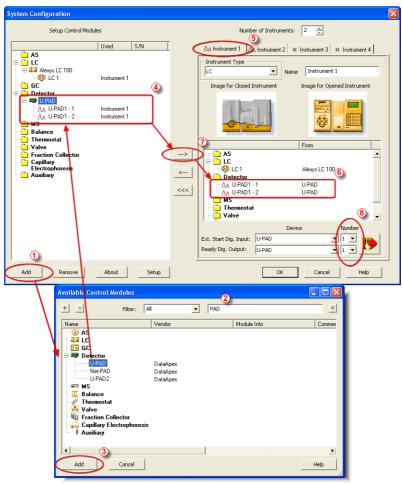


Fig. 8: System Configuration

- Start the Clarity station by clicking the icon on the desktop.
- Invoke the *System Configuration* dialog accessible from the *Clarity* window using the *System Configuration...* command.
- Press the Add button (1) (see Fig. 8 on pg. 9.) to invoke the Available Control Modules dialog.

- You can specify the searching filter ② to simplify the finding of the driver.
- Select the U-PAD and press the Add 3 button.

The DataApex U-PAD1 Setup dialog will appear.

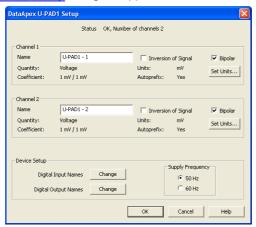


Fig. 9: DataApex U-PAD1 Setup

Enter the detector names in Name fields for individual channels, set signal units.

Note: A detailed description of this dialog can be found in the chapter "DataApex U-PAD1 Setup" on pg. 13.

• Press the OK button.

The **U-PAD** will appear in the Setup Control Modules list 4 of the System Configuration dialog.

- Drag and drop the U-PAD icon from the Setup Control Modules list 4 on the left side of the System Configuration dialog to the desired Instrument 5 tab on the right side 6 (or use the button 7 to do so).
- Set the Ext. Start Dig. Input and Ready Dig. Output numbers ⁽³⁾ for your acquisition card according to the wires being used for synchronization.

Note: The configuration dialog of the **U-PAD** device (<u>DataApex U-PAD1 Setup</u>) can be displayed any time by double-clicking on its icon or using the *Setup* button.

3.4.1 Measuring on multiple instruments

When measuring on multiple instruments using only a single **U-PAD** to gather data, rather than dragging the entire **U-PAD** icon from the *Setup Control Modules list* ①, drag the individual detector signals (**FID**, **UV-VIS** in this case) to separate Instruments ② (for more details see **Fig. 10** on pg. **11**.). The Instrument to which the particular signal is connected is listed to the right ③ of the detector name in the *Setup Control Modules* section.

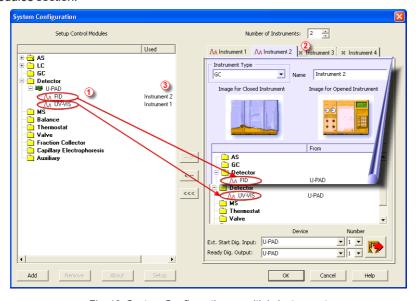


Fig. 10: System Configuration - multiple instruments

Note: Each independent Clarity must have separate Ext. Start Dig. Input and Ready Dig.

Output pins configured according to the actual wiring.

3.4.2 Using multiple U-PAD converters

When using multiple **U-PAD**s, after plugging the device into a USB port, the device will be detected automatically. After that, please, repeat the **U-PAD** configuration in **Clarity** described in the chapter **"Clarity Configuration"** on pg. 9.

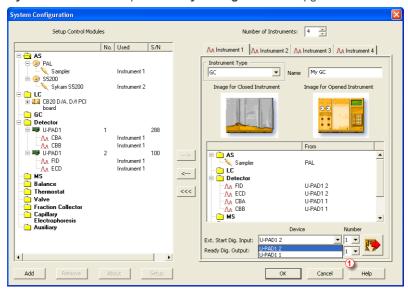


Fig. 11: System Configuration - multiple U-PADs

Individual converters can be assigned to multiple Instruments and combined together arbitrarily. In the example described in the **Fig. 11** on pg. **12**., all signals from **U-PAD1 1**, as well as all signals from **U-PAD1 2** are assigned to the *Instrument 1*.

While there are several cards assigned to the same Instrument, it is necessary to select which card will be used for *Ext. Start Dig. Input* and *Ready Dig. Output* signals ①.

All added A/D converters (e.g. **U-PAD**) must have their drivers properly installed. There must be correct number of items in the Windows OS *Device Manager*, otherwise you will not be able to add it in the *System Configuration* dialog and configure its detectors to Instruments.

Note: A maximum of **four U-PAD** converters can be configured and used in **Clarity** simultaneously.

Note:

A larger number of Instruments can be set than the amount you have purchased. You will not be able to measure on the surplus Instruments (indicated by a blue symbol of the curve on the tab), but you may use them e.g. for "offline" preparation of methods

4 Using the U-PAD

There are generally two places for setting the parameters of the **U-PAD A/D** Converter in Clarity:

- The <u>DataApex U-PAD1 Setup</u> dialog for setting the parameters that are set with respect to the type of connected signal.
- The Method Setup Acquisition tab for setting the parameters related to the type of analysis.

4.1 DataApex U-PAD1 Setup √ Full Version

The *DataApex U-PAD1 Setup* dialog sets the parameters of the **U-PAD** converter and since it depends on the type of connected signal, it is not going to be changed often.

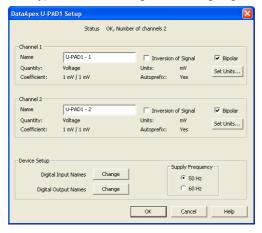


Fig. 12: DataApex U-PAD1 Setup

Number of Channels

Indicates the number of data acquisition channels.

Channel 1 (2)

For each channel of the **U-PAD A/D Converter** the name of the signal can be edited in the *Name* ① field and the *Set Units...* ② button can be used to change other signal parameters. Effects of these changes can be seen in other fields of the respective *Channel 1 (2) section.*

Name

This field will be used to describe the signal in the *Chromatogram* window.

Inversion of Signal

Inverts the polarity of the signal from the detector. This function can also be used for easy correction of inversely connected wires.

Bipolar

Both positive and negative voltages will be measured. Unchecking the *Bipolar* option will improve the resolution of the signal by 1 bit. On the other hand, it will also restrict the measuring of negative signals (i.e. below zero).

Set Units...

Invokes the *Detector Units* dialog that can be used for converting the voltage measured by **U-PAD** to the units of the detector. Such units will then be shown throughout **Clarity** instead of default mV. Setting is specific for each **U-PAD** channel (signal).



Fig. 13: Detector Units

- Quantity Field for entering a name of the y-axis in the graph of the measured signal.
- Units Field for entering the unit symbol, along with the prefix.
- Coefficient Allows to enter the coefficient between the selected units and mV (default).
- Autoprefix Allows to enter the unit prefix. In graphs, units will be scaled automatically, while in the tables, units with the prefix entered will be used.

Device Setup

This section allows to modify the naming of digital inputs and outputs and change the Mains Supply Frequency Suppression settings.

Digital Input Names

Allows to change the names of digital inputs.



Fig. 14: Digital Input Names

Digital Output Names

Allows to change the names of digital outputs. These names will be used in the *Digital Output Control* dialog in the chapter "Digital Inputs and Outputs" on pg. 19.



Fig. 15: Digital Output Names

Supply Frequency

Allows to change the Mains Supply Frequency Suppression. It is used for suppressing the interference on the specified frequency level. This setting influences the sampling frequency as well.

Note:

In most parts of Americas, the 60 Hz Mains frequency is usual, while the rest of the world tends to use 50 Hz Mains frequency (pre-set as default). The 50 Hz frequency is common in countries using 230V power supply, while the 60 Hz is connected to the 117V power supply.

4.2 Method Setup

Parameters of the *Method Setup - Acquisition* and *Method Setup - Measurement* dialogs depend on the type of analysis (measuring conditions), so they can be specific for various types of analyses. This is the reason for them to be a part of the method.

Caution: The parameters can't be changed during analysis.

4.2.1 Method Setup - Acquisition

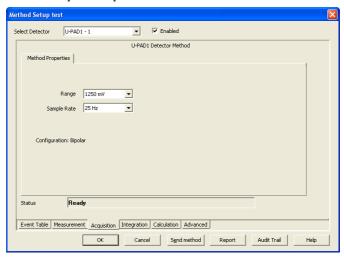


Fig. 16: Method Setup - Acquisition (U-PAD)

Select Detector

Selects the detector whose parameters are set. If there is only one detector (channel) configured on the **Clarity** Instrument, this control is not accessible.

Range

Selects the input voltage range (in mV). The selected range corresponds to the maximum output voltage of the detector that the A/D converter can still process. In case the detector unit conversion is set in Detector Units dialog, the range will be shown in detector units according to the conversion, while the corresponding original range in mV will be shown next to the field.

Available ranges for **U-PAD** converter are 156, 1 250 and 12 585 mV.

Sample Rate

Depending on *Supply Frequency* filter selected in the *System Configuration* – DataApex U-PAD1 Setup dialog, see the chapter "DataApex U-PAD1 Setup" on pg. 13. The available sampling rates are 6.25, 12.5, 25, 50 and 100 samples per second (for 50 Hz supply frequency filter) or 7.5, 15, 30, 60 and 120 samples per second (for 60 Hz supply frequency filter).

Higher sampling rate allows for the measurement of narrower peaks, but it also means a larger amount of data which affects the size of the resulting chromatogram and the speed of its processing.

Sample rate sufficient for successful peak detection is about 20 samples per narrowest peak.

A higher sampling rate also increases the amount of noise.

Note: Range and Sampling Rate are set individually for each detector.

4.2.2 Method Setup - Measurement

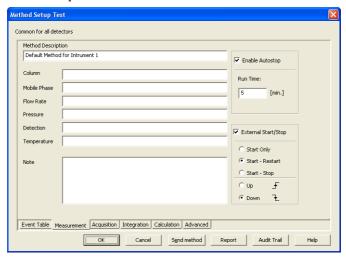


Fig. 17: Method Setup - Measurement

External Start/Stop

Enables control from an external signal.

Note:

The input used for an external start from a chromatograph can be set in the System Configuration dialog, see Fig. 8 on pg. 9.

Start Only

The external signal only starts an analysis. Triggering the input will not influence the acquisition in progress.

Start - Restart

The external signal terminates the analysis in progress and at the same time starts a new one. The first signal starts the analysis; the second signal terminates it and at the same time starts a new one - the so-called continuous measurement.

Start - Stop

The external signal starts and stops an analysis. The first signal starts the analysis; the second signal stops it, etc.

Down

The station reacts to changes in voltage in the relevant controlling input from high (> 3 V) to low (< 0.7 V), or to contacts closed in the relay.

Uр

The station reacts to changes in voltage in the relevant controlling input from low (< 0.7 V) to high (> 3 V), or to contacts opened in the relay.

4.3 Digital Inputs and Outputs

The **U-PAD A/D Converter** is equipped with two **digital TTL outputs**, which are also designed as relay contacts. It also has two **digital inputs** IN1 and IN2 equipped with optocouplers.



Fig. 18: Digital outputs of U-PAD1

These outputs can be controlled from the following places in Clarity:

• Standard assignment of external Ext. Start Dig. Input and Ready Dig. Output can be changed in the System Configuration dialog - 1 in the Fig. 8 on pg. 9.

Note: The assigned input and output is then automatically used in the ACTIVE sequence for synchronization with the autosampler.

- Digital outputs of U-PAD1 dialog.
- Device Monitor of the U-PAD1.
- In the Method Setup Event Table dialog.

4.4 Device Monitor

The *Device Monitor* window can be invoked by the *Monitor - Device Monitor* command from the *Instrument* window or using the **Device Monitor** icon. It displays the states of digital inputs and enables the user to monitor and change the states of digital outputs.



Fig. 19: Device Monitor - U-PAD1

Input no.

Lists the numbers of individual inputs.

Current State

Indicates the current status of the particular input. Green light corresponds to logical one (*HIGH*), light switched off (gray icon) to logical zero (*LOW*).

Descriptions

Shows the description indicating the meaning of individual input. The names of all digital inputs can be set in the DataApex U-PAD1 Setup dialog.

Output no.

Lists the serial numbers of individual outputs.

Current State

Indicates the current status of the particular output which also works as relays. Green light corresponds to logical one (*HIGH*), light switched off (gray icon) to logical zero (*LOW*). The state of any output can be changed by clicking on the *Current State* button of the particular output.

Descriptions

Shows the description indicating the meaning of individual output. The names of all digital outputs can be set in the DataApex U-PAD1 Setup dialog.

5 Troubleshooting

If you will not find your answers here, use the www.dataapex.com website where the Support menu will navigate you to frequently asked questions (FAQ), **Clarity** email conference archive or contact to **DataApex** helpdesk.

5.1 Locate your problem

When troubles occur, the fastest way to find a solution for it is to search for it in the following index via the **Dialog** (window), in which the problem occurred, **Error Messages** that appear or according to utilized **Hardware**. The name of the window/dialog is visible in its header.

Note:

Names of individual **Clarity Instruments** appear in the header instead of the common term "Instrument".

List of windows and dialogs

Dialogs	
Clarity	pg. 32 .
Data Acquisition	pg. 34 .
Instrument	pg. 32 .
Method Setup	pg. 32 .
Sequence	pg. 32 .
Single Analysis	pg. 32 .
System Configuration	pg. 22 ., pg. 32 ., pg. 34 .

List of Error messages

Ziot oi Ziroi moodagoo			
Error Messages			
Cannot establish communication with DataApex U-PAD	pg. 22 .		
Cannot load device driver	pg. 22 .		
Disabled (in the status line)	pg. 32 .		
Error occurred during setup	pg. 22 .		
Simulated (in Data Acquisition)	pg. 34 .		

Note:

To find solutions for error messages not listed above, check the **Clarity Getting Started** Guide or visit our web pages www.dataapex.com.

5.2 Problems with U-PAD

Most problems with **U-PAD** driver are generally indicated by following symptoms:

- U-PAD cannot be added to the Setup Control Modules in the System Configuration dialog, displaying "Cannot load device driver" error message followed by "Error occurred during setup".
- Claritywill issue an error message "Cannot establish communication with DataApex U-PAD" when opening the *Instrument* window.
- Double-clicking on the device will not invoke the <u>DataApex_U-PAD1_Setup</u> dialog.
- The Data Acquisition icon in the Instrument window is not active (this problem may also have another cause. For more details see the chapter Data Acquisition - non-functional on pg. 32.

Solution:

When one of these problems occur, first check if the **U-PAD** is correctly plugged in the **USB** port. If it is, check the **U-PAD** driver (see the chapter "How to check the **U-PAD** driver" on pg. 22.) and then proceed according to the suggested solutions

5.2.1 How to check the U-PAD driver

Windows XP/2000

In the lower left corner of the screen, in the *Start* menu, select *Settings - Control Panel* and the *System* icon, and under the *Hardware* tab click on the *Device Manager* button.

Windows 7/Vista

In the lower left corner of the screen, in the *Start* menu, type "Device Manager" in the search field and click on the found item.

Windows 8.1

Press the *Windows* key on the keyboard and start typing *Device Manager*, in the search results, click on the desired item to display the *Device Manager* window.

In the desired *Device Manager* window look up the resulting list for item "Chromatography Devices - U-PAD v. 1.3 (DataApex Ltd.)".

In the following image, all drivers (except the **U-PAD 2.0** driver) are installed correctly.

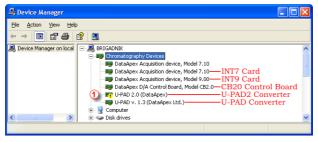


Fig. 20: Device Manager in Windows XP

Error status of the driver may be one of the following:

If the "Chromatography Device – U-PAD v. 1.3 (DataApex Ltd.)" - item does not appear, then the driver has not been installed.

Solution:

Reinstall the driver manually as described in the chapter "Installation and reinstallation in Windows XP/Vista/7/8.1" on pg. 24.

The driver is installed, but there is a yellow exclamation mark (1) in the Fig. 20 on pg. 22.) or a question mark over its icon, or it is not behaving as expected. The driver is probably installed incorrectly.

Solution:

Reinstall the driver manually as described in the chapter "Installation and reinstallation in Windows XP/Vista/7/8.1" on pg. 24.

The item does appear, but is displayed with a cross symbol.

Solution:

Double-click the item to invoke the *General* tab and activate the driver by clicking on the *Enable this device* button.

The item does appear and there are no error symbols around it. A driver for a different type of Windows operating system was possibly installed thereby replacing the correct driver.

Solution:

Uninstall the incorrect driver and install the correct one as described in the chapter "Installation and reinstallation in Windows XP/Vista/7/8.1" on pg. 24.

5.3 Manual installation

5.3.1 Installation and reinstallation in Windows XP/Vista/7/8.1

- Connect the U-PAD with a cable to a USB port on the computer.
- During start-up, Windows should automatically recognize the new Plug and Play device and start the Add Hardware Wizard.

Note: If the wizard does not occur, use the Start - Control Panel - Add Hardware to invoke installation.



Fig. 21: Step 1 of Hardware Installation Wizard

Click the Next button.



Fig. 22: Step 2 of Hardware Installation Wizard

• Select No, not this time and click Next.

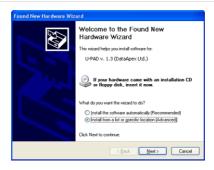


Fig. 23: Step 3 of Hardware Installation Wizard

• Select the Install from a list or specific location (Advanced) option and click Next.



Fig. 24: Step 4 of Hardware Installation Wizard

• Select Don't search. I Will choose the driver to install, and click the Next button.

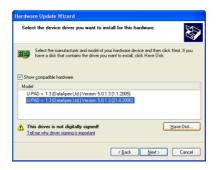


Fig. 25: Step 5 of Hardware Installation Wizard

• To be sure that **Windows** install the correct driver, click the *Have Disk...* button.



Fig. 26: Step 6 of Hardware Installation Wizard

Click the Browse... button and in the displayed Locate File dialog browse to the
main folder of the Clarity Station and HW_DRIVERS\UPAD subfolder
(C:\CLARITY\HW_DRIVERS\UPAD by default). Select the CSWUPAD.INF file
and click the Open button. Now you are back in the Install From Disk dialog
displaying the correct path to the CSWUPAD.INF file. Click the OK button.

Note: The driver is also located on the Clarity installation DVD-ROM in the HW_DRIVERS\UPAD subfolder.

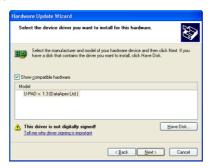


Fig. 27: Step 7 of Hardware Installation Wizard

 Now the Hardware Update Wizard shows the correct driver to be installed in Windows. Click the Next button and after completing the installation click the Finish button in the following dialog.

5.3.1.1 Reinstallation of drivers using the System Restore Point

Use the **System Restore Point** in **MS Windows XP/Vista/7** to uninstall the incorrect drivers

- Go to Start Programs Accessories System Tools System Restore in Windows XP or press the Windows key on your keyboard and type "System Restore" in the search field and click on the found item in Windows Vista/7/8.1.
- In the System Restore dialog use the Restore my computer to an earlier time in Windows XP or Choose a different restore point in Vista/7 option and press the Next button.
- In the next dialog select a restore point from a time before the incorrect driver was installed.

Caution: Any software or hardware that was installed after the moment of the selected restore point will be removed from your system.

- Reinstall Clarity and then restart the PC.
- Windows XP should now correctly identify and install the drivers.

5.3.2 Reinstallation of drivers in Windows 2000

Note: Windows **2000** is supported until **Clarity** version **3.0** only.

After installing **Clarity** and the **U-PAD** device, the *Device Manager* either does not display the *Chromatography Devices* – *U-PAD* item at all or it displays a yellow exclamation mark next to the item that indicates that the device is not working properly.



Fig. 28: Device Manager in Windows 2000

If the "U-PAD v. 1.3 (DataApex Ltd.)" item is missing, Windows has automatically installed incorrect driver.

If you are able to locate which device was mistakenly installed instead of the **U-PAD** driver, you can open its *Properties* and *Update this driver*. In the manual installation process, select the correct drivers analogically to the description in the chapter "Installing the correct driver manually" on pg. 29.

If the correct driver cannot be determined, the last installed driver will have to be removed.

Caution: This procedure can be used safely only if you did not install any other hardware since the installation of the converter.

5.3.2.1 Deleting incorrect *.INF files in the system

Note: Be sure to have the **Clarity**software already installed.

Open the C:\WINNT\INF (C:\WINDOWS\INF) folder ① in Windows Explorer.

Note: You may have to allow for the hidden system files to be displayed. In menu select Tools – Folder Options - View tab then set the Hidden Files and Folders option to Show Hidden Files and Folders.

- Sort the files according to Modified column ② and locate the last modified *.PNF file ③.
- Also locate the corresponding *.INF file 3 and delete both of them.



Fig. 29: Windows Explorer

 Continue to install the correct driver manually (described in the chapter "Installing the correct driver manually" on pg. 29.) without restarting the PC.

5.3.2.2 Installing the correct driver manually

The following procedure described step by step can be applied under any circumstances.

- Invoke the Start Control Panel Add/Remove Hardware command.
- In the first dialog press the Next button.



Fig. 30: Step 1 of Add/Remove Hardware Wizard

• Select Add/Troubleshoot a device and click the Next button.



Fig. 31: Step 2 of Add/Remove Hardware Wizard

• In the **Devices** list select the *Add a new device* and click the *Next* button.



Fig. 32: Step 3 of Add/Remove Hardware Wizard

• Select No. I want to select the hardware from a list and click the Next button.



Fig. 33: Step 4 of Add/Remove Hardware Wizard

 In the Hardware types list select the Other devices option and press the Next button.



Fig. 34: Step 5 of Add/Remove Hardware Wizard

Set the path to the correct drivers. These drivers can be found in the directory
where Clarity is installed (C:\CLARITY\BIN by default), in the HW_
DRIVERS\UPAD subfolder. Then click the OK button.

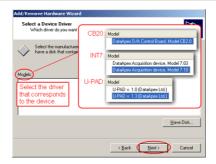


Fig. 35: Step 6 of Add/Remove Hardware Wizard

- In the Models list select the U-PAD v. 1.3 (DataApex Ltd.) and click the Next button.
- In the remaining dialogs press the Next button.
- · Restart the PC.
- Re-check the driver in the *Device Manager* as described in the chapter "Reinstallation of drivers in Windows 2000" on pg. 27.

5.4 Data Acquisition - non-functional

Gray icon with the heading DISABLED and non-functional Monitor - Data Acquisition command.



Fig. 36: Data Acquisition disabled

Other manifestations of this error are also: <u>Method Setup - Acquisition</u> tab missing, *Method - Acquisition* command non-functional, *Run*, *Stop*, *Abort* and other similar commands non-functional in the *Single Analysis*, *Sequence* or *Device Monitor* windows. Possible causes are:

Description:

a) You are using **Clarity Offline** or a **Clarity DEMO** version, which does not enable the measurement of chromatograms.



Fig. 37: Clarity Offline and Clarity DEMO

Solution:

Check whether there is a blue line with the title **OFFLINE** displayed in the main *Clarity* window under the symbols of the Instruments, or the title **DEMO** in the *Clarity* window header.

In case of Clarity DEMO, use the Windows Start menu to locate the Clarity DEMO group in the *All Programs* section and select the *Remove Clarity* item there.

Then, use the correct **Clarity** installation DVD to install the full version.

In case of the **Clarity Offline** station, remove the HW key with the **Offline** license and insert key with the **Clarity** full licence.

Description: b) Detector is not allocated to the Clarity Instrument:



Fig. 38: Detector not allocated to the Instrument

Open the *System Configuration* dialog from the *Clarity* window using the *System - Configuration...* command and check the tab of the corresponding Instrument - **Instrument X**. If it has no allocated detectors, add them.

Solution:

In the left-hand list Setup Control Modules select the correct detector connected to the A/D card you are using and drag it to the corresponding instrument on the right.

If your **A/D card** is not in the left-hand list *Setup Control Modules*, add it using the *Add* button and repeat the previous step.

Note:

More information on the *System Configuration* dialog can be found in the chapter "Clarity Configuration" on pq. 9.

Description:

c) You have a licence purchased for data collection from a smaller number of Instruments:

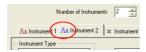


Fig. 39: Small number of Instruments purchased

Open the *System Configuration* dialog from the *Clarity* window using the *System-Configuration...* command and check the tab of the corresponding Instrument - **Instrument X**. Blue symbol of the curve in the tab header indicates that Instrument is not usable for measuring.

Solution:

Check your serial number S/N for example using the command *Help - About...* from the main *Clarity* window.

5.5 Data Acquisition - Simulated

The title "Simulated" is displayed in the Data Acquisition window.



Fig. 40: Data Acquisition Simulated

Description: The corresponding Instrument only displays the simulated curve (from the file

 ${\it CHANNX.DTA file}) in the \textit{\it Data Acquisition} \ window, or in other words \ has \ allocated$

a **DEMO detector** on itself.

Solution: Open the System Configuration dialog from the Clarity window using the System -

Configuration... command and check the tab of the corresponding Instrument - Instrument X. If it only has the detector signals from the DataApex DEMO detector, it is necessary to reconfigure it. You can find more information how to

add a detector in the chapter "Clarity Configuration" on pg. 9...

Make sure that you have installed a full version of **Clarity**. If this is not the case, inscription *DEMO* is in the header of the main window. If so, install the full version and enter the user code you have been provided. Also make sure that the detector

is properly configured.

6 Tables and specifications

6.1 Description of the INT7 connector (Male)

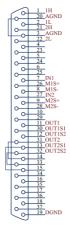


Fig. 41: Connector pins

Tab. 3: Description of the connector pins:

Pin	Description
1H	+ Input of the 1st channel
1L	- Input of the 1st channel
AGND	Ground of the analog channels
IN1	Digital input (of the 1st channel). Used as an External Start (marker) of the 1st channel
OUT1	1st Digital output Digital outputs are controlled by Event Table, ACTIVE sequence and Digital Output Control UPAD dialog.
DGND	Ground of the digital inputs and outputs.
M1S+	+ pin of the 1st Digital input. Level is <i>LOW</i> when current flows between M1S+ and M1S
M1S-	- pin of the 1st Digital input.
OUT1S1 OUT1S2	Digital outputs - 1st relay contact. Contact is closed when output is at a <i>LOW</i> level.

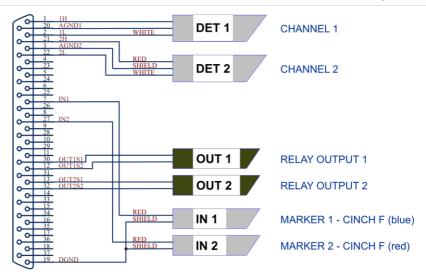


Fig. 42: Standard U-PAD cable for two detectors

6.2 Parameters of digital inputs and outputs

Digital Outputs

The board contains two digital TTL outputs (OUT1 and OUT2) with maximum permissible current of 20 mA at both levels. Both digital outputs are also designed as relay contacts (OUT1S1/OUT1S2 and OUT2S1/OUT2S2).

The maximum permissible current is 100~mA, the maximum-switched voltage is 100~V, and the contact resistance is $150~m\Omega$. The relay is on-state (closed) at level LOW of a relevant digital output.

Digital Inputs

The **U-PAD** A/D converter contains two digital TTL inputs IN1, IN2. The inputs are also equipped with an optocoupler (M1S+/M1S- and M2S+/M2S-). The excitation current is 5 *mA* and corresponds to input level *LOW*. The LED of the isolator is directly connected to the connector.

Caution: A resistor must be inserted (e.g. in the connector proper) to restrict the diode current to 10 mA max.

To suppress contact bouncing the digital inputs are equipped with monostable toggles that prolong any change affecting the input to at least 200 ms (to ensure that the station is able to detect it). A pulse with duration of at least 1 ms is necessary to excite the toggles.

6.3 U-PAD - CE Conformity Declaration

CE CONFORMITY DECLARATION

We, DataApex Ltd.

Podohradska 1704/1

CZ - 155 00 Prague 5

declare under our sole responsibility that

the product:

chromatography integrator

type:

U-PAD

to which this declaration relates is in conformity with the following standards

EN 61326-1:98+Z1:99+Z2:02

following the provision of Directive:

89/336/EEC including amendments

Complementary information: Certificate EZÚ no.: 1030724 dated 23.4.2003

The last two digits of the year in which the CE marking was affixed: 03

Place of issue: Prague Manufacturer representative: L. Siska



Date of issue: 6.5.2003 Position: hardware production manager

Fig. 43: Conformity Declaration

6.4 Technical data

Tab. 4: Technical data of the U-PAD device:

Parameter	Value
Converter type:	24-bit δΣ (delta – sigma)
Analog inputs:	differential

Parameter	Value			
Number of channels:	2, isolated from USB (inputs share common ground)			
Input range:	unipolar and bipolar, each channel independently 156, 1 250 or 12 585 mV			
Input impedance:	1 ΜΩ			
Sample rate:	6.25, 12.5, 25, 50, 100 samples per second for 50 Hz mains frequency filter 7.5, 15, 30, 60, 120 samples per second for 60 Hz mains frequency filter			
Non-linearity:	< 0.003 %			
Digital inputs:	2 TTL (5 V), can be also controlled using the built-in optocouplers ((M1S+/M1S- and M2S+/M2S-, 5 mA current)			
Digital outputs:	$2\ TTL\ (5\ V,\ maximum\ current\ load\ 20\ mA),\ also\ as$ relay contacts (switching contacts rating: max. switching current 0.5 ADC, max. contact resistance $150\ m\Omega,\ max.\ voltage\ 100\ V)$			
Dimensions:	100 x 55 x 16 mm			
Weight:	110 g			
Power supply:	from the PC via the USB cable			

Tab. 5: Effective resolution for various acquisition speeds and input ranges:

	12 585 mV	1 250 mV	156 mV
6.25 Hz	22 bit	22 bit	20 bit
12.5 Hz	21 bit	21.5 bit	20 bit
50 Hz	20.5 bit	21 bit	19 bit
100 Hz	20 bit	20 bit	18.5 bit

6.5 Disposal and Recycling



Electrical and electronic equipment must **not be** disposed of as unsorted municipal waste.

It should be collected separately to ensure proper treatment and recycling. Please return used equipment to designated collection points.