



SHIMADZU LC-10/20 PUMP

Clarity Control Module

ENG

Code/Rev.: M091/90F

Date: 2026-05-22

Phone: +420 251 013 400

clarity@dataapex.com

www.dataapex.com

DataApex Ltd.
Petřilkova 2583/13
158 00 Prague 5
Czech Republic

Clarity[®], DataApex[®] and ▲[®] are trademarks of DataApex Ltd. Microsoft[®] and Windows[™] are trademarks of Microsoft Corporation.
DataApex reserves the right to make changes to manuals without prior notice. Updated manuals can be downloaded from www.dataapex.com.

Author: DM

Contents

1 Shimadzu LC-10/20 Pump	1
2 Requirements	2
3 Installation Procedure	3
3.1 Shimadzu LC-10/20 Pump communication	3
3.2 Clarity Configuration	5
4 Using the control module	7
4.1 Method Setup - LC Gradient	7
4.1.1 Gradient Options	10
4.2 Method Setup - Advanced	11
4.3 Device Monitor	12
4.3.1 LC Control Manual Flow	14
4.4 Shimadzu LC-10/20 Pump Setup	15
5 Report Setup	17

To facilitate the orientation in the **Shimadzu LC-10/20 Pump** manual and **Clarity** chromatography station, different fonts are used throughout the manual. Meanings of these fonts are:

Open File (italics) describes the commands and names of fields in **Clarity**, parameters that can be entered into them or a window or dialog name.

WORK1 (capitals) indicates the name of the file and/or directory.

ACTIVE (capital italics) marks the state of the station or its part.

Chromatogram (blue underlined) marks clickable links referring to related chapters.

The bold text is sometimes also used for important parts of the text and the name of the **Clarity** station. Moreover, some sections are written in format other than normal text. These sections are formatted as follows:

Note: Notifies the reader of relevant information.

Caution: Warns the user of possibly dangerous or very important information.

Marks the problem statement or trouble question.

Description: Presents more detailed information on the problem, describes its causes, etc.

Solution: Marks the response to the question, presents a procedure how to remove it.

1 Shimadzu LC-10/20 Pump

This manual describes the setting of the **Shimadzu LC-10/20 Pump**, which may be used to create high-pressure or low-pressure gradient. The control module enables direct control of the **LC-10ADvp**, **LC-10ATvp** and **LC-20A** (whole series, **LC-10** emulation mode) over serial line.

Caution: The **Shimadzu LC-10/20 Pump** control module can not be combined with **Shimadzu LC-10/20 System** control module on a single PC where Microsoft .NET Framework is installed.

The **Shimadzu LC-10/20 Pump** control module can control only a single pump on a single PC, which means that it is not possible to control more than one (of above mentioned) pump on a single PC through this control module.



Fig. 1: Shimadzu LC-10ADvp Pump

Direct control means that the pump can be completely controlled from the **Clarity** environment. Instrument method controlling the analysis conditions will be saved in the measured chromatograms.

2 Requirements

- **Clarity** Installation with LC Control module (p/n A24).
- Free USB port in the PC.
- Opt-USB convertor (p/n OPT01).

Note: The **DataApex Opt-USB** converter conforms to the similar **Shimadzu** converter (Shimadzu p/n 22-35480-91).

- Pump firmware version 2.2 or greater.

Caution: Minimal required firmware version for **LC-10ADvp** is **5.26** and for **LC-10ATvp** is **5.27**.

3 Installation Procedure

3.1 Shimadzu LC-10/20 Pump communication

The **Shimadzu LC-10/20 Pump** is controlled by serial (RS232) communication. The communication is provided via the Optical to USB converter connected to the computer via USB A-A cable (p/n SK10) and to the pump via optical cable (p/n SK15). Both of these cables are part of the **DataApex Opt-USB** converter set (p/n OPT01), but can also be ordered separately.

Optical to USB converter

For communication an OPT-RS232 converter is needed. It can either be the **DataApex Opt-USB** converter (p/n OPT01) or **Shimadzu** adaptor (Shimadzu p/n 22-35480-91).

Pump Settings

Communication parameters should be set on the pump in the *SERVICE MODE* (see pump documentation) as follows:

Baud rate 2400, parity N, bits 8, stop bit 1.

SYS (Set System Parameter) has to be set to 5 for isocratic mode or to 6 for gradient mode from the pump keyboard.

.....
Caution: These values are not documented in the pump manual.

The other Pump system parameters should be set to:

Local=0 (Remote)

ADRS= any value (not used)

CBM LINK=0 (not used)

Shimadzu LC-20 settings

Shimadzu LC-20 pumps have to be set to the **LC-10** emulation mode in order to be able to work in **Clarity**. This is done from the pump's keyboard according to this procedure:

- Press the **VP** button on the pump's front keyboard until the *CALIBRATION* inscription is shown on the pump's display.
- Press the **func** button to enter the pump's configuration. The password has to be filled in.

.....
Note: The default password is 00000.

- Press the **func** button repeatedly until the *OP MODE* inscription is shown on the pump's display.
- Enter the desired number (1) to switch the pump to the **SCL-10Avp** emulation mode and press the **enter** button.
- Switch the pump off and on again. The new communication parameters should be saved and used.

The procedure is in detail described in the **LC-20** hardware manual provided by the **Shimadzu** company.

3.2 Clarity Configuration

- In the *System Configuration* dialog press the *Add* button (① on Fig 5) to invoke the *Available Control Modules* dialog.

Fig. 2: Available Control Modules

- Select the **Shimadzu LC-10/20 Pump** in the **LC** section and press the *Add* (② on Fig 5) button.
- The [Shimadzu LC-10/20 Pump Setup](#) dialog will appear.

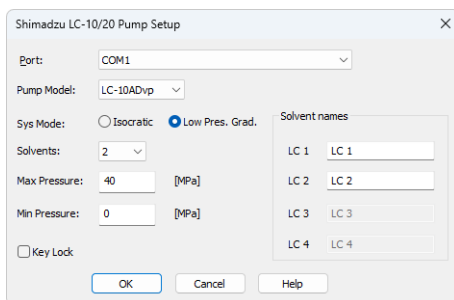


Fig. 3: Shimadzu LC-10/20 Pump Setup

- Select the correct communication *Port* and fill the other necessary parameters in this dialog.

Note: Other fields and parameters of the [Shimadzu LC-10/20 Pump Setup](#) dialog are described later in the manual (for more details see the chapter **Shimadzu LC-10/20 Pump Setup** on pg. 15)..

- The **Shimadzu LC-10/20 Pump** item (③ with the number of solvents set in the [Shimadzu LC-10/20 Pump Setup](#) dialog will appear in the *Setup Control Modules* list of the *System Configuration* dialog.
- Change the *Instrument Type* (⑤ a) on the desired *Instrument* tab (⑤ b) to LC.
- Drag the LC item from the *Setup Control Modules* list on the left to the *Instrument* tab on the right (⑥), or use the (→) button (⑦).

Note: The **Shimadzu LC-10/20 Pump** can be used for the low-pressure gradient with the use of switching valves as the pump can handle up to 4 solvents.

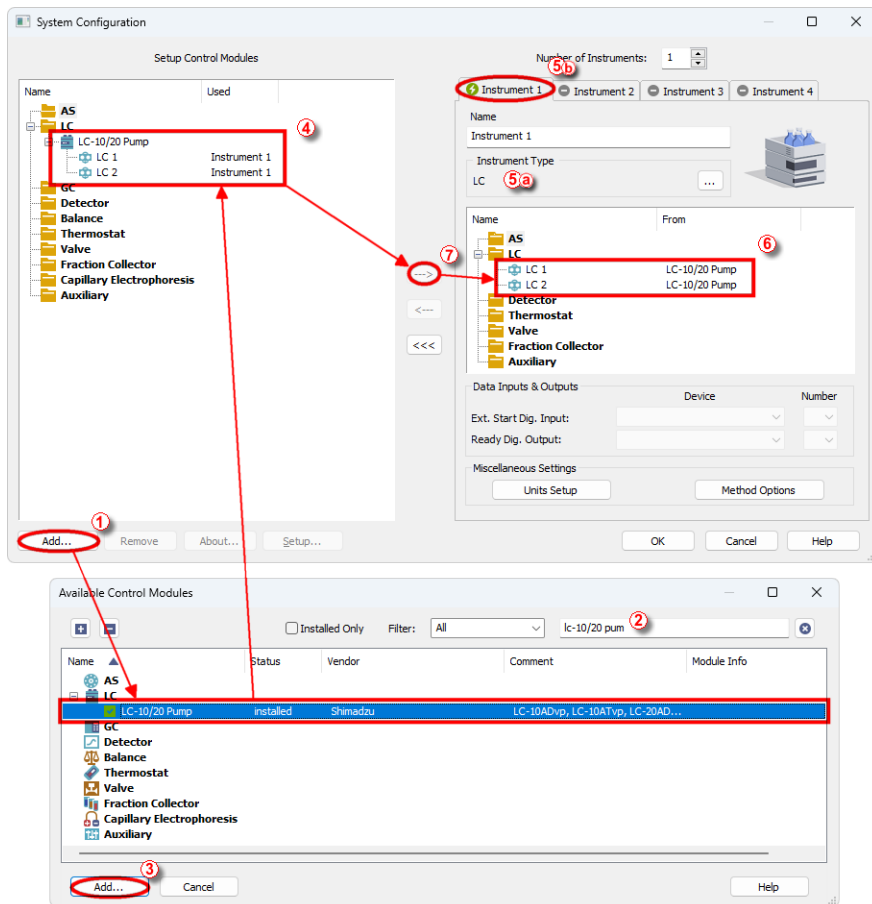


Fig. 4: System Configuration

4 Using the control module

New [LC Gradient](#) tab appears in the *Method Setup* dialog, enabling the setting of the LC gradient program.

4.1 Method Setup - LC Gradient

The *Method Setup - LC Gradient* dialog serves for setting up the LC instrument method.

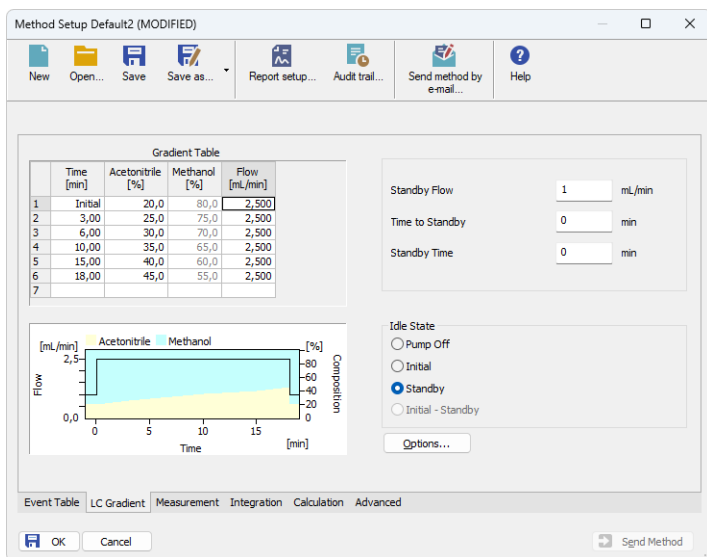


Fig. 5: Method Setup - LC Gradient

Gradient Table

A table for setting the composition of the mobile phase and the overall flow rate as a function of time. Operation is analogous to that of spreadsheets (Excel, Quatro Pro, etc.). To prepare the cell to receive values, click it by the left mouse button; the cell will highlight by dots. A cell that fails to highlight is not available for editing.

Time [min.]

Sets the time at which the ratio of flow rates and the overall flow rate correspond to the values entered in the corresponding row. (These values vary continuously from one time to the next in a manner ensuring that the conditions specified in the next row are satisfied).

XXX1 (..4) [%]

Represents the percentage of a component. The designation **XXX1-4** is in fact replaced by the name of the component (items *Solvent 1 - 4* in the [Gradient Options](#) dialog). Should you enter a component value such that the sum of all values exceeds 100 %, the percentage in the last column is automatically

adjusted; if the percentage of the last compound is already zero, the value of the currently entered component is adjusted instead. The flow rate of a compound is calculated by multiplying the overall flow rate (indicated in the *Flow* column) by the corresponding percentage divided by 100.

Flow [ml/min]

Indicates the overall flow rate through the column. The entered value applies to the time specified in the corresponding row.

Graph

The graph depicts the percentage of components as a function of time together with the overall flow rate. Data are taken over from the **Gradient Table**. Changes effected in this table are immediately reflected in the graph. Legend in the header of the graph indicates the assignment of colors to individual components. The assignment is fixed and individual components are displayed in the graph from bottom to top. The flow rate is displayed as a black line.

The graph has two vertical axes: the axis on the left refers to the mixing ratio, the one on the right to the overall flow rate.

Parameters

Standby Flow

Sets the overall flow rate through the column in the *STANDBY* state reached after the last row of the table has been performed and the time period defined in the *Time to Standby* field has passed. The duration of this state is defined by the *Standby Time* item. The ratio of individual components in the respective *STANDBY* and *IDLE* states is given by the first row of the **Gradient Table** (the *Initial* row).

Time to Standby [min]

Indicates the time during which the flow rate and mobile phase composition changes continuously between the last values entered in the table and the values defined by *Standby Flow* field and the *Initial* row mobile phase composition.

This time is included in the analysis time (the Instrument is in the *CONTROL* state). In case when the *Time to Standby* is zero, there is step change from flow and components percentage specified on the last row of gradient table to that specified for *STANDBY* state.

Standby Time [min]

The time during which the flow rate is maintained at *Standby Flow*. This time is included in the analysis time (the Instrument is in the *CONTROL* state).

Initial - Standby

The **Shimadzu LC-10/20 Pump** control module doesn't support this option.

Idle State

An item specifying the overall flow rate through the column outside the instrument method. The following options are possible:

Pump Off

The flow rates of all components are zero.

Caution: Be careful as this setting may damage the column in some cases.

Initial

The flow rate is defined by the first row of the gradient table (the *Initial* row).

Standby

The flow rate is the same as in the *STANDBY* mode and, accordingly, corresponds to the value entered in *Standby Flow* field.

The *IDLE* state comes into effect each time an Instrument is opened, at the end or after abortion of an analysis by the *Abort* command, and is also maintained after the **Clarity** program is shut down.

The mixing ratio of individual components in both the *IDLE* and *STANDBY* states is given by the first row of the **Gradient Table** (the *Initial* row).

Note: There is a step change in the flow and components percentage from the values specified for the *STANDBY* state to those specified for the *IDLE* state if the *Idle State* field is not set to *Standby*.

4.1.1 Gradient Options

Invoke the *Options...* button in the [Method Setup - LC Gradient](#) dialog to open the *Gradient Options* dialog. This dialog allows to set the custom name for particular solvents, to switch whether they are used or not in the gradient and to set the warning levels for pressure to prevent the damage to hardware.

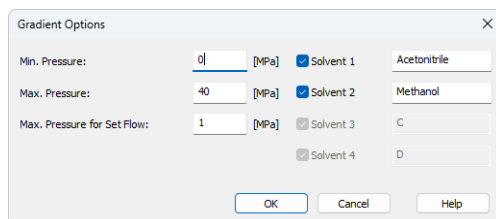


Fig. 6: Gradient Options

Min. Pressure

Sets the minimum pressure for the given pump. In case of a low-pressure condition, a 1-minute timeout is applied to allow the pump to re-pressurize before shutdown is triggered to prevent solvent leakage. The pressure limit can be also set in [System Configuration - Shimadzu LC-10/20 Pump Setup](#) dialog. The minimum pressure defined in the method cannot exceed the limit specified in *System Configuration*.

Max. Pressure

Sets the maximum pressure for the given pump. When pressure reaches the set value, the pump will shut down to prevent the damage to the pump when the column is blocked. The pressure limit can be also set in [System Configuration - Shimadzu LC-10/20 Pump Setup](#) dialog. The maximum pressure defined in the method cannot exceed the limit specified in *System Configuration*.

Max. Pressure for Set Flow

Sets the maximum pressure to be used in the *Set Flow* dialog in the [Device Monitor](#). Initial value is 1 MPa.

Solvent 1 (.4)

It is possible to enable/disable particular solvent, as well as to set custom name to it.

4.2 Method Setup - Advanced

Method Setup Default2 (MODIFIED)

New Open... Save Save as... Report setup... Audit trail... Send method by e-mail... Help

Common for all detectors

Subtraction
 Chromatogram [None]
 Matching No Change
 Set... None

Column Calculations
 Unretained Time 0 [min]
 Column Length 50 [mm]
 Statistical Moments
 From Width at 50%

	Auxiliary Signal	Store
1	LC Pressure LC 1	<input checked="" type="checkbox"/>
2	LC Flow LC 2	<input checked="" type="checkbox"/>

User Variables
 Variable 1
 Name MethodUserVar1
 Value 0
 Variable 2
 Name MethodUserVar2
 Value 0
 Variable 3
 Name MethodUserVar3
 Value 0


Event Table LC Gradient Measurement Acquisition Integration Calculation **Advanced**

OK Cancel Send Method

Fig. 7: Method Setup - Advanced

Shimadzu LC-10/20 Pump provides the pump flow and pressure auxiliary signals for use in **Clarity**. To save the auxiliary signals into each chromatogram measured according to the given method, check the particular checkbox in the lower section of the Method Setup - Advanced dialog.

4.3 Device Monitor

The pump status dialog can be invoked by the *Monitor - Device Monitor* command from the *Instrument* window or using the Device Monitor  icon. It displays the actual flows of particular solvents, as well as the total flow, the total pressure and the analysis time. In the individual section of the **Shimadzu LC-10/20 Pump** the status of the pump may be seen.

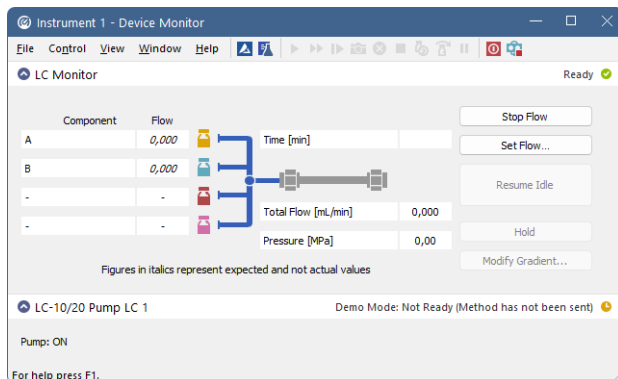


Fig. 8: LC Monitor

Stop Flow

The pumps can be stopped from this window using the *Stop Flow* button. This action will stop the pump only, the analysis run will continue and must be stopped or aborted separately by the *Stop/Abort* button in the toolbar.

Set Flow...

Sets the desired total flow and solvent ratios in the opened *Set Flow* dialog. *Max. Pressure* allows to set a different pressure than set in the "Gradient Options" on page 10 of a given method.

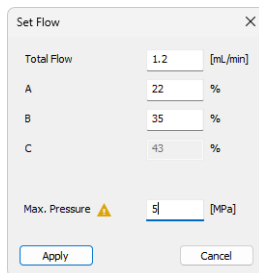


Fig. 9: Set Flow

Resume Idle

Returns the pumps to *IDLE* state as defined in the appropriate field on the [LC Gradient](#) tab of the *Method Setup* dialog.

Hold/Resume

Clicking on the *Hold* button will keep the current gradient conditions until *Resume* is clicked.

Modify Gradient...

Opens the [LC Control Manual Flow](#) dialog allowing to set custom flow and mobile phase composition, disregarding the **Gradient Table** set in the method. The command is only available during the analysis run.

4.3.1 LC Control Manual Flow

The *LC Control Manual Flow* dialog accessible through using the *Modify Gradient ...* button from [Device Monitor](#) window allows the user to set a custom mobile phase composition and flow while the analysis is running. It resembles the [Method Setup - LC Gradient](#) tab in functionality.

The screenshot shows the 'LC Control Manual Flow' dialog box. It features a 'Gradient Table' with the following data:

	Time [min]	Acetonitrile [%]	Methanol [%]	Ethanol [%]	Flow [mL/min]
1	Initial	80,0	20,0	0,0	2,500
2	3,00	80,0	20,0	0,0	2,500
3	6,00	60,0	40,0	0,0	2,500
4	10,00	60,0	40,0	0,0	2,500
5	15,00	20,0	80,0	0,0	2,500
6	18,00	20,0	80,0	0,0	2,500

Below the table is a graph showing 'Flow [mL/min]' on the left y-axis (0,0 to 2,5) and 'Composition [%]' on the right y-axis (0 to 80) against 'Time [min]' on the x-axis (0 to 15). The graph shows a constant flow of 2.5 mL/min. The composition starts at 80% Acetonitrile and 20% Methanol, then at 6 minutes, it changes to 60% Acetonitrile and 40% Methanol, and at 15 minutes, it changes to 20% Acetonitrile and 80% Methanol. The legend indicates Acetonitrile (yellow), Methanol (cyan), and Ethanol (pink).

Control settings on the right include:

- Standby Flow: 1 mL/min
- Time to Standby: 0 min
- Standby Time: 0 min
- Idle State:
 - Pump Off
 - Initial
 - Standby
 - Initial - Standby

Buttons for 'OK' and 'Cancel' are located at the bottom right.

Fig. 10: LC Control Manual Flow dialog

The *LC Control Manual Flow* dialog is only available during the analysis run. When it is invoked and the **OK** button is pressed, the original **Gradient Table** from the acquisition method is discarded and replaced by the **Gradient Table** from the *LC Control Manual Flow* dialog. Any such operation is recorded in the audit trail of the measured chromatogram.

Caution: After the analysis run which used manual flow changes ends, the original method is automatically sent to all controlled devices to make sure **Clarity** station returns to the original method.

4.4 Shimadzu LC-10/20 Pump Setup

Shimadzu LC-10/20 Pump Setup dialog (accessible through the *System Configuration* dialog) is used to set the connection to the pump, select its type and set some other basic parameters.

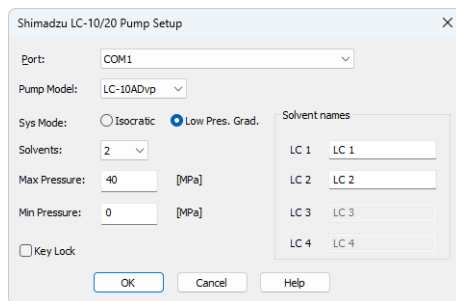


Fig. 11: Shimadzu LC-10/20 Pump Setup

Port

Sets the COM port to which the **Shimadzu LC-10/20 Pump** pump is connected.

Pump Model

Allows to select from several types of Shimadzu pumps. **LC-10ADvp** and **LC-10ATvp** models are supported at the moment.

Note: If any **LC-20** type pump should be connected, it is necessary to set the **LC-10 emulation mode** on the pump (for more details see the chapter **Installation Procedure** on pg. 3)..

Sys Mode

Sets the pump operation mode. *Isocratic* and *Low Pres. Grad.* options are available.

Note: The pump hardware must be set accordingly (for more details see the chapter **Installation Procedure** on pg. 3)..

Solvents

Sets the number of solvents when the *Low Pres. Grad.* mode is used.

Max Pressure

Sets the maximum pressure for the given pump. When pressure reaches the set value, the pump will shut down to prevent the damage to the pump when the column is blocked. The pressure limit can be also set in [Method Setup - LC Gradient - Gradient Options](#) dialog. The maximum pressure defined in the method cannot exceed the limit specified in *System Configuration*.

Min Pressure

Sets the minimum pressure for the given pump. In case of a low-pressure condition, a 1-minute timeout is applied to allow the pump to re-pressurize before shutdown is triggered to prevent solvent leakage. The pressure limit can be also set in [Method](#)

[Setup - LC Gradient - Gradient Options](#) dialog. The minimum pressure defined in the method cannot exceed the limit specified in *System Configuration*.

Key Lock

Serves for making the **Shimadzu LC-10/20 Pump** keyboard inactive.

5 Report Setup

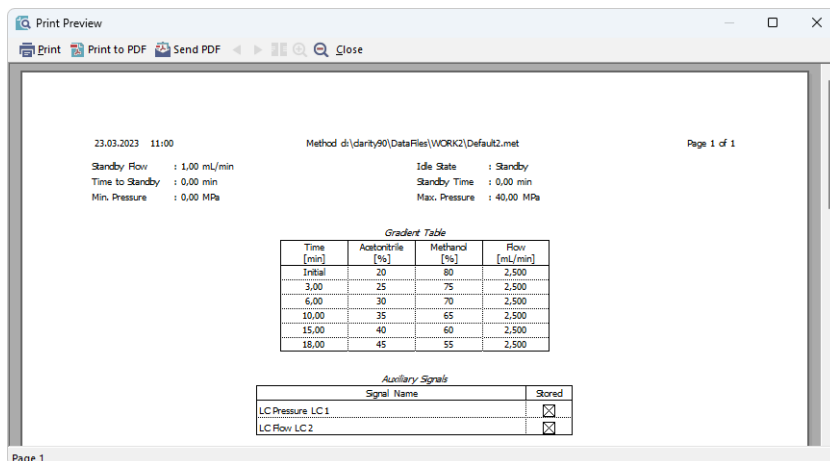


Fig. 12: Shimadzu LC-10/20 Pump report preview

All general LC settings (such as the **Gradient Table** from the [Method Setup - LC Gradient](#) tab) are reported as a part of the data displayed by the use of *Instrument Control* checkbox of the *Report Setup - Method* dialog.