



CTC PAL3 SYSTEM

Clarity Control Module

ENG

Code/Rev.: M224/80F
Date: 2/16/2021

Phone: +420 251 013 400
Fax: +420 251 013 401
clarity@dataapex.com
www.dataapex.com

DataApex Ltd.
Petrzilkova 2583/13
158 00 Prague 5
The Czech Republic

Clarity[®], DataApex[®] and ▲[®] are trademarks of DataApex Ltd. Microsoft[®] and Windows[™] are trademarks of Microsoft Corporation.

DataApex reserves the right to make changes to manuals without prior notice. Updated manuals can be downloaded from www.dataapex.com.

Author: LD

Contents

1 CTC PAL3 System	1
2 Requirements	2
2.1 Software requirements	2
2.2 Hardware requirements	2
2.3 Limitations	2
3 Installation procedure	3
3.1 Installing Correct Version of Agilent ICF	3
3.2 Network connections	4
3.3 Clarity Configuration	6
3.4 Installation Qualification of Agilent ICF	10
4 Using CTC PAL3 System	12
4.1 Device Monitor	12
4.2 Method Setup	14
4.3 Method Setup - Advanced	17
5 Troubleshooting	18
5.1 Specific Problems	19

To facilitate the orientation in the **CTC PAL3 System** manual and **Clarity** chromatography station, different fonts are used throughout the manual. Meanings of these fonts are:

Instrument (blue text) marks the name of the window to which the text refers.

Open File (italics) describes the commands and names of fields in **Clarity**, parameters that can be entered into them or a window or dialog name (when you already are in the topic describing the window).

WORK1 (capitals) indicates the name of the file and/or directory.

ACTIVE (capital italics) marks the state of the station or its part.

The bold text is sometimes also used for important parts of the text and the name of the **Clarity** station. Moreover, some sections are written in format other than normal text. These sections are formatted as follows:

Note: Notifies the reader of relevant information.

Caution: Warns the user of possibly dangerous or very important information.

■ Marks the problem statement or trouble question.

Description: Presents more detailed information on the problem, describes its causes, etc.

Solution: Marks the response to the question, presents a procedure how to remove it.

1 CTC PAL3 System

This manual describes the use of the **PAL RTC**, **PAL LSI** and **PAL RSI** autosamplers with the **Clarity** software ver. **8.0 and later**. **Clarity** comes with ICF version **A.02.04 Revised**.



Fig 1: CTC PAL3 System

The control module enables direct control of the instrument over Local Area Network (LAN). Direct control means that the autosampler can be completely controlled from the Clarity environment. The Instrument method controlling the analysis conditions will be saved in the measured chromatograms.

2 Requirements

2.1 Software requirements

Clarity (p/n C50) Installation USB with **AS control** module (p/n A26).

Agilent ICF and **CTC PAL3 System** require **Microsoft .NET version 4.0** for correct installation. This version is already installed on majority of PCs. Nonetheless you will be notified during the installation if your PC is missing any version of **Microsoft .NET** - then follow the instructions there. Required operating systems:

- **Windows 7 SP1 (32/64 bit)**
- **Windows 8.1 (32/64 bit)**
- **Windows 10 (32/64 bit)**

For complete list of .NET requirements, see the **.NET Framework System Requirements** on Microsoft web page.

Caution Before installing **Clarity**, make sure that your **Windows** are updated to the latest version.

2.2 Hardware requirements

- Version of firmware must be compatible with the **Agilent ICF** installed with **Clarity**.
- LAN interface installed on PC is required.

2.3 Limitations

- Help is not fully functional in all windows.
- Demo mode is not supported.
- Injection overlapping for headspace sampler is not implemented yet.

3 Installation procedure

CTC PAL3 System is not part of the **Clarity Typical** installation. To install it, select the *Full* installation or *Custom* installation with ICF option selected.

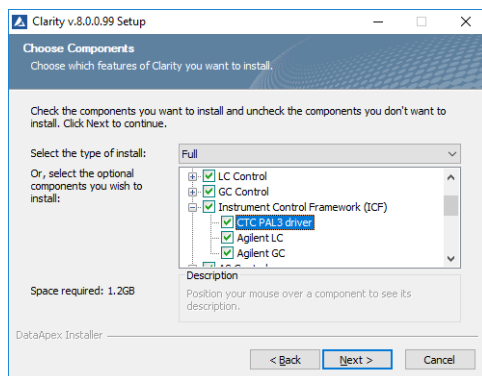


Fig 2: Full installation of Clarity

3.1 Installing Correct Version of Agilent ICF

Clarity expects a specific version of Agilent ICF. Because other programs may also be using Agilent ICF, it is possible they've installed a different version than is supported by Clarity. In that case in the installation you will be prompted to reinstall it in order to install correct version. Going forward with this step is crucial for correct functionality of Clarity and Agilent ICF. This situation may also occur during Clarity update when the new version contains updated ICF version.

Note: This reinstallation of Agilent ICF may cause that other programs using it, may not function properly.

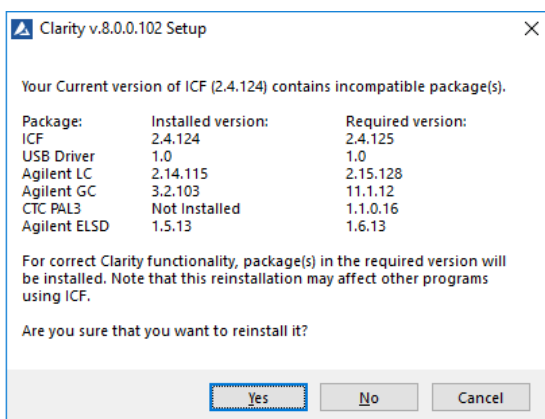


Fig 3: Installing correct version of Agilent ICF

3.2 Network connections

The instruments supported by **Agilent ICF** has to be connected to a site network by LAN. It is recommended to attach the hardware directly to the PC avoiding hubs, switches etc. When using a switch or a hub, multiple hardware can be connected to one PC. Always contact your local LAN administrator who can make the appropriate settings.

Caution: **Cross LAN** cable is primarily used for the direct connection of the instrument and the PC. This cable can also be used for the connection of the device to the switch or network socket, but with older switches, the **straight LAN** cable might be necessary.

LAN Settings

PC: LAN card, TCP/IP protocol.

Both PC and supported CTC hardware should be configured on the same IP range.

Tab 1: Recommended IP settings:

	CTC Hardware	LAN card
IP	192.168.111.235	192.168.0.1
Port	64000	- - -
Subnet Mask	255.255.255.0	255.255.255.0
Default Gateway	192.168.0.1	192.168.0.1
Accept From IP	0.0.0.0	- - -
Accept From Subnet Mask	0.0.0.0	- - -

Firewall

Ensure that the firewall does not block communication from the **CTC** hardware.

Internet connection

Be aware that many Internet connections are provided using the LAN card. If this is the case a separate LAN card must be configured for each Internet and connected **CTC** hardware.

3.3 Clarity Configuration

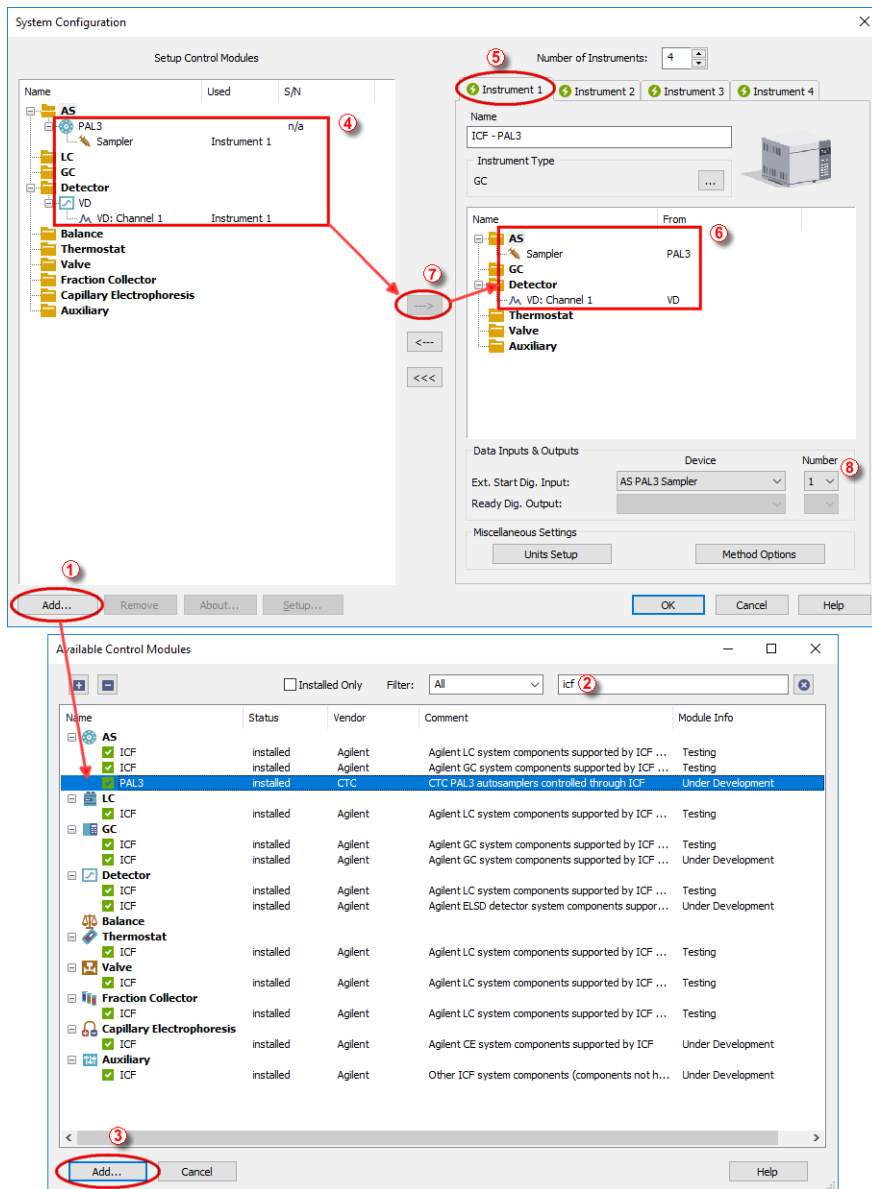



Fig 4: System Configuration

Caution: Before you start **Clarity**, ensure there is not any other application active and controlling the instruments.

- Start the **Clarity** station by clicking on the  icon on the desktop.
- Invoke the **System Configuration** dialog accessible from the **Clarity** window using the *System - Configuration...* command.
- Press the **Add** button ① (See **Fig 4** on pg 6.) to invoke the **Available Control Modules** dialog.
- You can specify the searching filter ② to simplify the finding of the driver.
- Select the **PAL3** item and press the **Add** ③ button.

The PAL3 Setup dialog will appear.

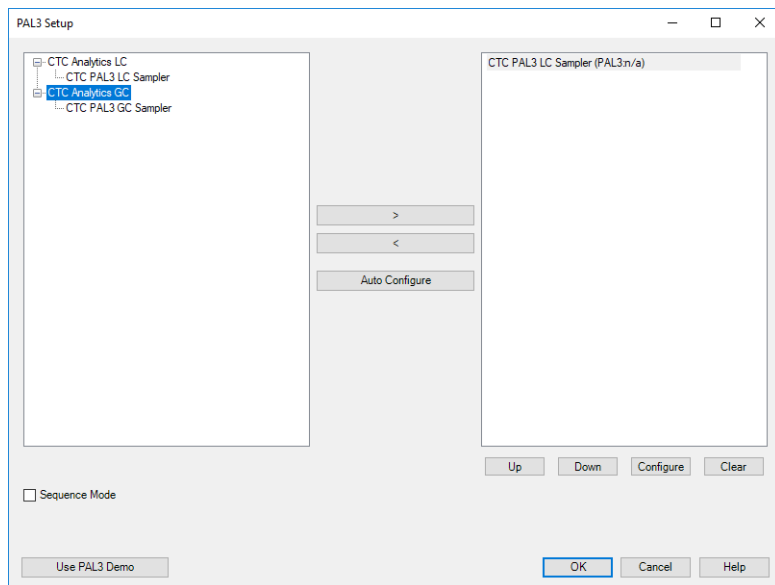


Fig 5: PAL3 Setup

Note: Press the *F1* key to display the **CTC PAL3 System** help with detailed description of the dialog.

- Click on the *Auto Configure* button in the **PAL3 Setup** dialog and the **Automatic configuration parameters** dialog is displayed for entering the connection parameters.

Note: *Auto Configure* button is active only when CTC Analytics LC or CTC Analytics GC item is selected.

Note: The *Sequence Mode* option can optimize timing for maximum of 50 measurements. In case there are more measurements in sequence the optimization of sequence will be splitted into groups containing maximum 50 measurements.

Note: CTC PAL3 GC Sampler should be used to run GC PAL3 Scripts. CTC PAL3 LC Sampler should be used to run LC PAL3 Scripts.

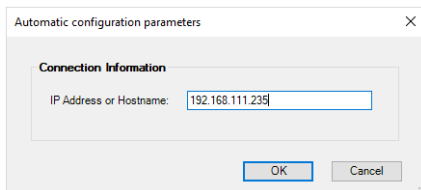


Fig 6: Automatic configuration parameters dialog

- Fill in the *IP address* or *Hostname* and click *OK* button.

For manual configuration click on *Configure* button. Sampler's *Configure* dialog will appear.

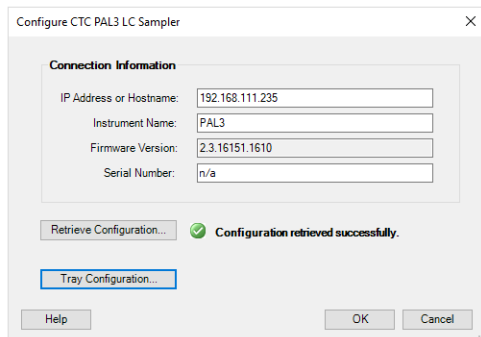


Fig 7: Automatic configuration parameters dialog



Note: The user has to fill the serial number manually. The serial number is not stored in the PAL firmware.

Note: Configuration can be retrieve using the *Retrieve Configuration* button. If the device is offline the answer is: *Retreive configuration failed. No response from Server...* If the device is already used by other user the answer is: *Retreive configuration failed. The role is already taken.*

- Fill in required parameters and click *OK* button.
- Then click *OK* button in the lower part of the [PAL3 Setup](#) dialog.

The **CTC PAL3 System** will appear in the *Setup Control Modules* list ④ of the [System Configuration](#) dialog.

Caution: When the *PAL3 Setup* dialog is open for extended time, communication between **Clarity** and the AS may be lost, in such case an error message will be displayed. After accepting the error message it is necessary to set the configuration from very beginning, starting with clicking the *Add* button ① (See **Fig 4** on pg 6.) in the **System Configuration** dialog.

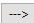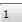- Drag and drop the **PAL3** icon from the *Setup Control Modules* list ④ on the left side of the **System Configuration** dialog to the desired *Instrument* ⑤ tab on the right side ⑥ (or use the  button ⑦ to do so).
- Set the *Ext. Start Dig. Input* ⑧ to "1" -  ⑧ for ICF Sampler to assure correctly synchronized start of the analysis with the **PAL3** sampler's injection.

Caution: If you change the configuration of the **CTC PAL3 System** (for example remove and add another instruments), the communication parameters like *IP address* will be lost and it is necessary to set them again.

Note: The configuration dialog of the **CTC PAL3 System** can be displayed any time by double-clicking on its icon or using the *Setup* button.

3.4 Installation Qualification of Agilent ICF

Agilent ICF is an external program developed by Agilent and for that reason it must be validated using their utility. If you have installed Clarity with Agilent ICF, **the IQ is valid only if successful validation of ICF is attached.**

The validation of ICF can be performed directly from the [IQ Report](#).

Caution: If Clarity expects ICF installation then **IQ** also. And if for some reason the ICF installation is not found, the *Installation Qualification Test* is automatically **FAILED**. To resolve this situation, it is necessary to re-install **Agilent ICF** during Clarity installation.

Date	27.03.2018, 16:28
Serial number of application	000-00000+000/255
User Code	Serial number could not be found.
Version of application	Clarity version 8.0.0.67
Build date of application	25.03.2018, 23:37
Instruments	
Extensions	
Controls	
Certification file	C:\Clarity\Bin\iq.chk
Checksum of cert. file	
Date of cert. file	26.03.2018, 01:27
User	kana
System	Microsoft Windows 10 Professional version 10.0 (Build 16299)
Acquisition and hardware devices	Key Rockey 4ND 088-00888-38910/000 id:179B4ECF ICF CN10939118

Installation Qualification Test: Passed

⚠ Agilent ICF is present in your system, it is necessary to perform its validation separately, the IQ result is valid only if successful validation of ICF is attached. To perform validation [click here](#) ①

Files

File	Path	Version	Size	File date	Status
AdvionCMS.dll	C:\Clarity\Bin	5.0.20.1	385536	26.03.2018, 01:11	Passed
advioncommon.dll	C:\Clarity\Bin	5.0.20.1	16384	26.03.2018, 01:11	Passed
AdvionData.dll	C:\Clarity\Bin	5.0.20.1	321536	26.03.2018, 01:11	Passed
AdvionObjects.dll	C:\Clarity\Bin	5.0.20.1	626176	26.03.2018, 01:11	Passed
ASPEN.dll	C:\Clarity\Bin	1.2.4.0	185344	26.03.2018, 01:11	Passed
asp64.sys	c:\clarity\bin\hw_drivers\asp2	4.7.0.8	84832	26.03.2018, 01:11	Passed

Fig 8: IQ Report with ICF installation present

Press the link "click here" ① and after that click the *Run* button to generate reports. They will be automatically opened in the in your default internet browser. There is a total number of 5 reports.

Address field of the generated report displays the location of the actual report ②.

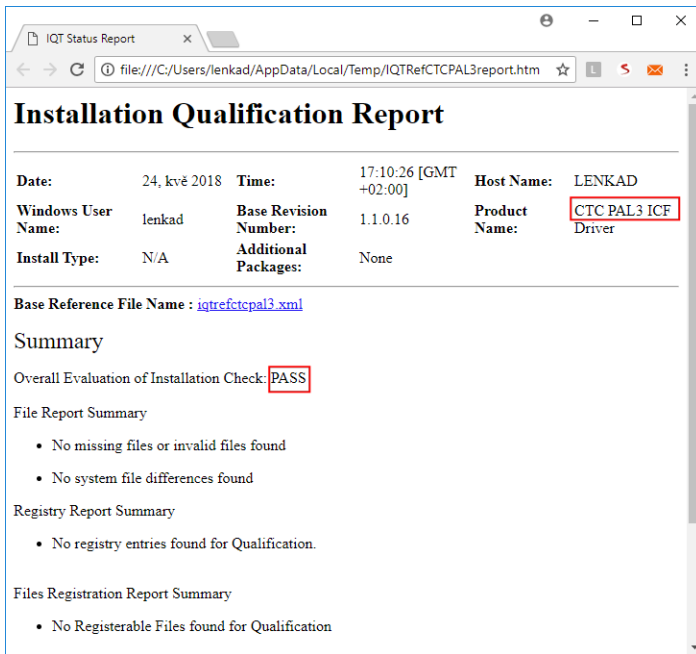


Fig 9: ICF Report - PASS


4 Using CTC PAL3 System

There are multiple places for setting the parameters of the **CTC PAL3 System** in the **Clarity**:

- the [Device Monitor](#)
- the [Method Setup](#)

Caution: Before opening the [Instrument](#) window with configured **ICF** devices, ensure there is not any other PC connected to AS. Otherwise there will be raised an error during the connection.

4.1 Device Monitor

Before using CTC PAL3 System, corresponding scripts must be loaded from the [Device Monitor](#) window using the Script Manager. The [Device Monitor](#) window can be invoked by the *Monitor - Device Monitor* command from the [Instrument](#) window or using the **Device Monitor**  icon. Then Script Manager can be invoked from **Device Monitor** by holding right mouse button over PAL3 schema for 2-3s.

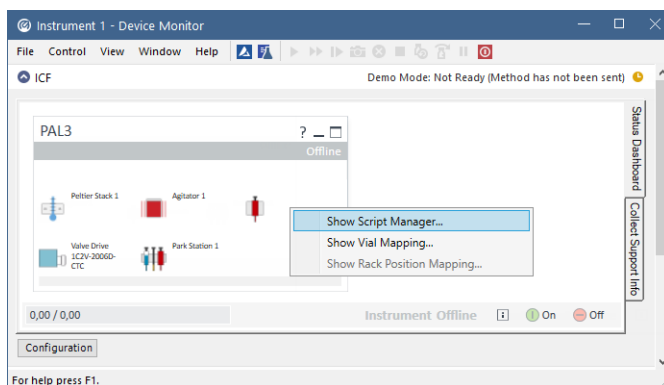


Fig 10: Device Monitor

Use *Import Scripts* button and select the available scripts from `Clarity\Bin\HW_DRIVERS\CTC_PAL3_Scripts`.

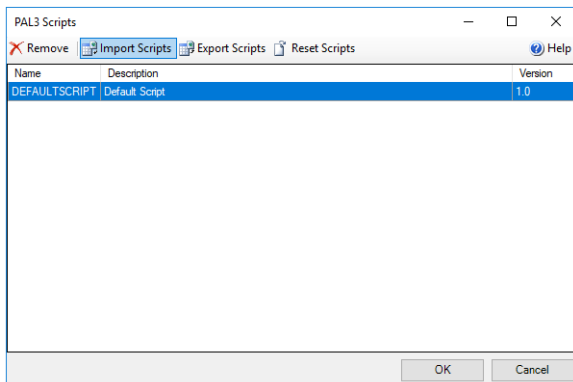


Fig 11: Script Manager

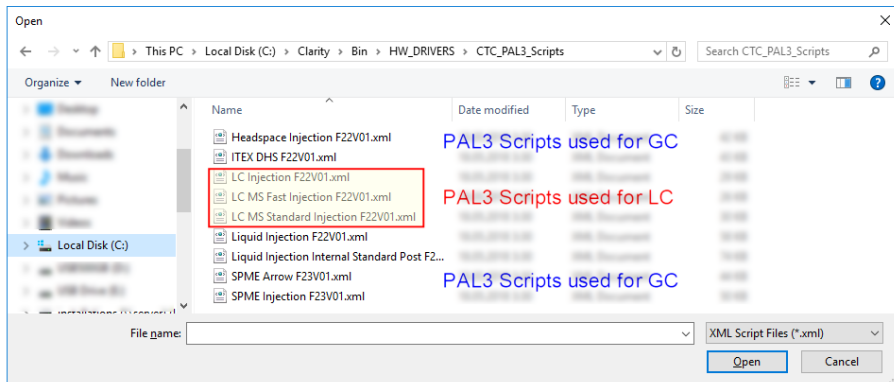


Fig 12: Script selection

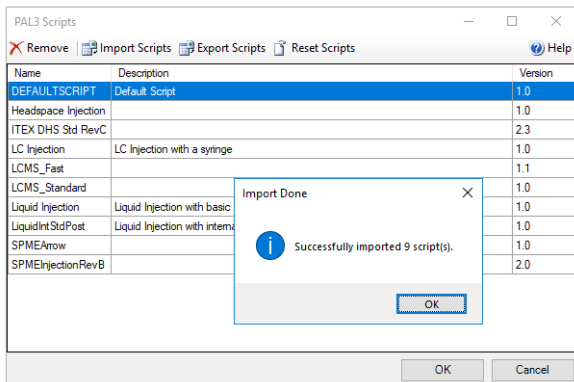


Fig 13: Import confirmation

Button *Configuration* on the bottom part of autosampler section in the *Device Monitor* allows to enter and modify configuration of autosampler directly from the *Device Monitor*. This feature is intended for quick updating of configuration when trays or other tool of the autosampler are changed. Modifications of configuration may cause that method might need an adaptation before its use in analysis. The *Configuration* feature is not accessible during running analysis or running *Sequence*.

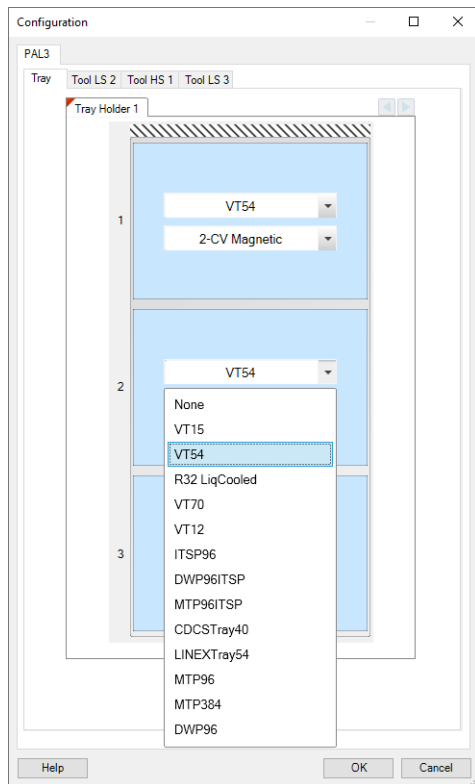


Fig 14: Configuration from Device Monitor

4.2 Method Setup

In **AS** tab you can select appropriate Method Script ① and then set all the parameters for AS control ②.

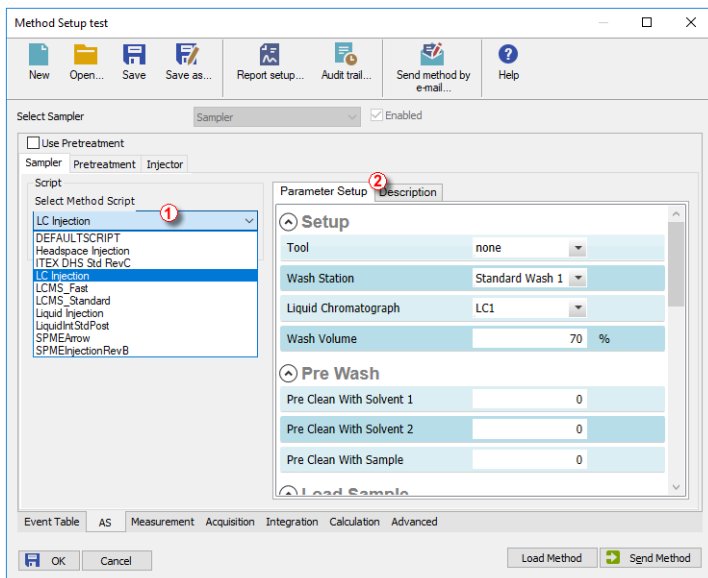


Fig 15: Method Setup - Parameter Setup

Description tab is mainly informational.

Pretreatment tab is not supported.

Injector tab is not editable, it displays the injector selected by the script.

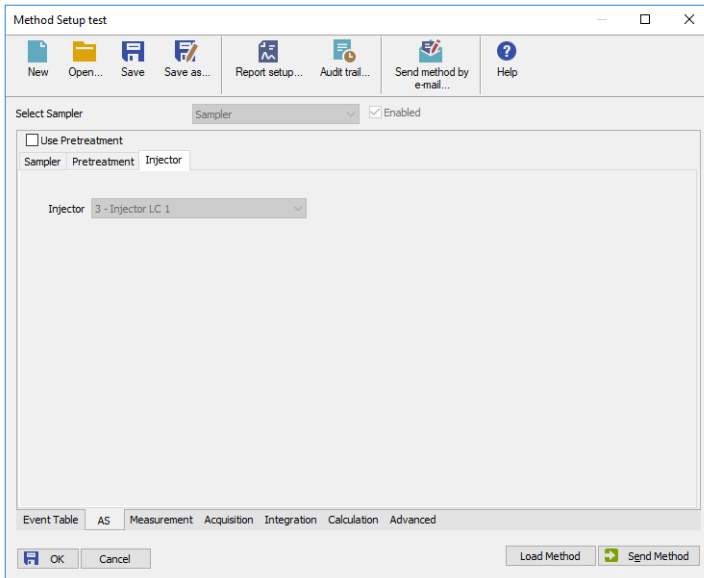


Fig 16: Method Setup - Injector

4.3 Method Setup - Advanced

The **Method Setup - Advanced** tab serves for setting the usage of auxiliary signals of the **CTC PAL3 System** autosampler.

Method Setup Default1

Common for all detectors

Subtraction
Chromatogram:
Matching:

Column Calculations
Unretained Time: [min]
Column Length: [mm]
 Statistical Moments
 From Width at 50%

	Auxiliary Signal	Store
1	Agitator 1 - PAL3	<input type="checkbox"/>
2	SyringeHeater - PAL3	<input type="checkbox"/>

User Variables
Variable 1
Name:
Value:
Variable 2
Name:
Value:
Variable 3
Name:
Value:

Event Table AS Measurement Acquisition Integration Calculation **Advanced**

Fig 17: Method Setup - Advanced

The list of available auxiliary signals is shown in the table in the lower part of the dialog. By checking the checkbox in the *Store* column for the particular row, the given auxiliary signal will be stored into the measured chromatogram.

Note: LC Injection valve auxiliary signal serves for service purposes only, under standard conditions it gives no data.

5 Troubleshooting

When the remedy for some problem cannot be discovered easily, the recording of communication between **Clarity** and **CTC PAL3 System** control module can significantly help the **DataApex** support to discover the cause of the problem.

The recording can be enabled by adding or amending the LOGGING.INI file in the **Clarity** installation directory (C:\CLARITY\CFG by default). The file can be edited in any text editor (e.g. Notepad). Following section should be edited:

```
[Log]
echo = ON
filename = log_%D.txt
reset = OFF
; Sections List:
AuditTrail = ON
BadTrace = ON
CommandLine = ON
Acquisit = ON
AgilentICF = ON
SST = OFF
Internet = OFF
FractionCollector = OFF
```

Note: %D (or %d) in the filename parameter means that the log will be created separately for each day. The *reset = OFF* parameter disables deleting the content of the log each time the station is started during the same day.

The created *.TXT files will greatly help in diagnosis of unrecognized errors and problems.

In case you cannot establish communication with Agilent ICF controlled instruments, please review the following issues:

Check the network connection using the Ping command


The problem in communication between **Clarity** and Agilent instruments may be caused by wrong network configuration, firewall preventing the connection, etc. Run the command line in Windows (for example by pressing the **Windows key** together with the **R** key, in the displayed **Run** window type *cmd* and press *Enter*).

In the command line type ping <ip-address-of-instrument> or <hostname> and press *Enter*. The *IP Address (hostname)* is the same you entered in the ICF;Setup dialog.

5.1 Specific Problems

Clarity can't be run and it displays "Agilent ICF is not installed correctly." message.

Cause: The cause of the problem is that the Agilent ICF has a different version than expected by Clarity. It can typically happen when other software also using Agilent ICF decides to reinstall it. Thus next time Clarity expects different version than is installed.

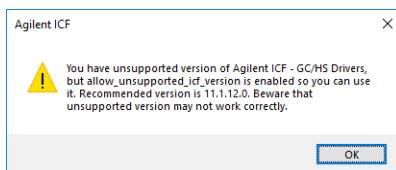

Solution: Solution is to reinstall Agilent ICF during Clarity installation. Follow steps described in the chapter "**Installation procedure**" on pg 3.

Analysis can't be run in headspace mode and Clarity displays "Voltage of PWMOut is not set to 34V" message.

Cause: In the Headspace mode, the settings of the PWMOut are expected to be set to 34V through the controller in order to be able to switch the valve. Default value after installation is 0V (for non-headspace samplers) in order not to power the valve which does not have to be present.

Solution: Solution is to switch the PAL3 using the controller through *Setup - PAL Modules - I/O Interfaces - PWMOut - Output Voltage* to 34V. Please note that in order to be able to set the value, the user must be switched to the "Extended User" access level (done by simultaneously pressing both buttons on the controller).