

Clarity2Go

Analysis monitoring for smartphones

Clarity2Go is a mobile companion to Clarity Chromatography Station.

It enables you to monitor selected parameters of running sequences or single analyses conveniently from your mobile. Its use is free for all users of Clarity.

The monitoring works only if both the smartphone and the PC with Clarity have access to the internet.

The whole solution consists of three parts:


- Clarity Software (Full, minimal version 4.0.4) communicating with the web server
- Clarity2Go web server run by DataApex (provided to Clarity2Go users free of charge)
- Clarity2Go application for smartphones and tablets

Clarity2Go application is currently available for iOS ([Apple Store](#)) and Android ([Google Play](#)).

How to use Clarity2Go

Enabling Clarity2Go in Clarity

- In Clarity use *System – Clarity2Go...* to invoke *Clarity2Go Configurations* dialog.
- In the dialog, check the Instruments that you want to monitor by Clarity2Go. In the example below, Instruments 1 and 2 will be monitored.
- It is possible to protect the monitoring by a password which can be set in *Advanced >>* options.

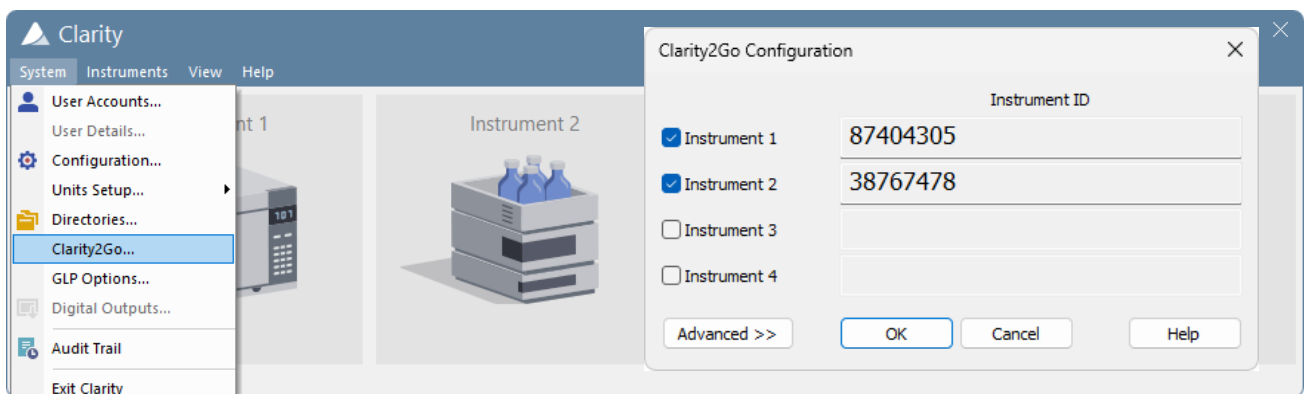

- Enter the *Instrument ID* (and the optional password) in Clarity2Go and you are ready to monitor the state of your instrument.

Clarity2Go on your smartphone and tablet

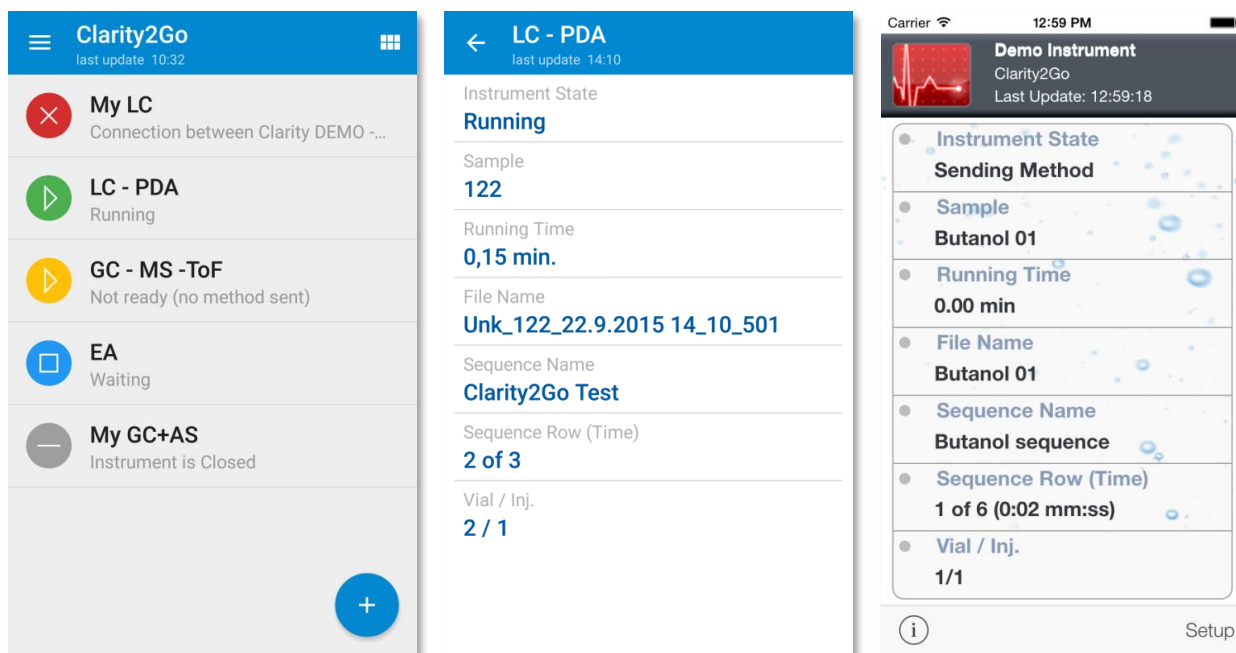


Figure 1. Clarity2Go on smartphones with Android (first two images) and with iOS (last image)

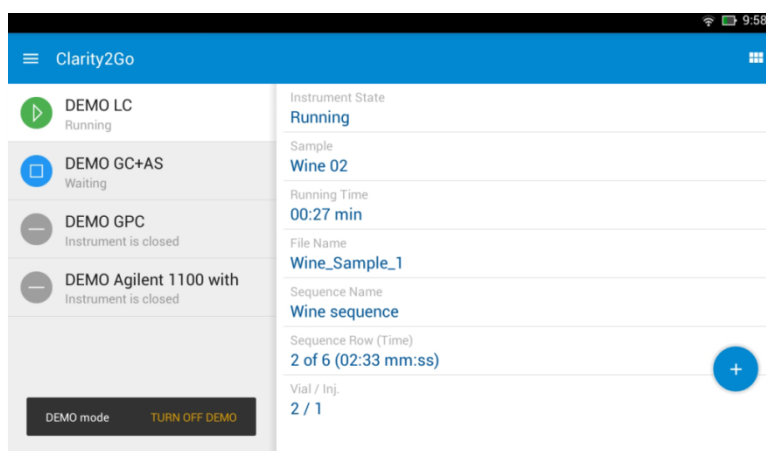


Figure 2. Clarity2Go on an Android tablet

The development of Clarity2Go is currently being put on hold; if you encounter any problems, please contact support@dataapex.com.