

## Multilingual Clarity (p/n C50) and Clarity Lite (p/n C40)

Clarity (Lite) Chromatography Software is supplied with multiple languages from version 5.0.


Available languages: **English, Spanish, French, Russian, Chinese** and **German**. German is freely available from Clarity (Lite) version 5.0.4.

Translation is limited to the main SW, Extensions and most of Control Modules. Control modules localization might differ for each language and DataApex will be adding individual localized drivers in coming Clarity versions. Help and documentation are in English and will be also localized continuously. All data files are independent on the language and can be used in any localization.


## How to get Multilingual Clarity (Lite)

### New users

New licenses are provided in the latest available version so users will automatically get the multilingual Clarity. You can set the preferred language at the start of the installation:

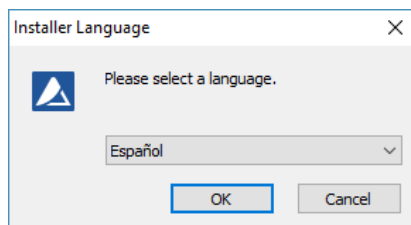

Or anytime later using the *Help – Languages* menu command:

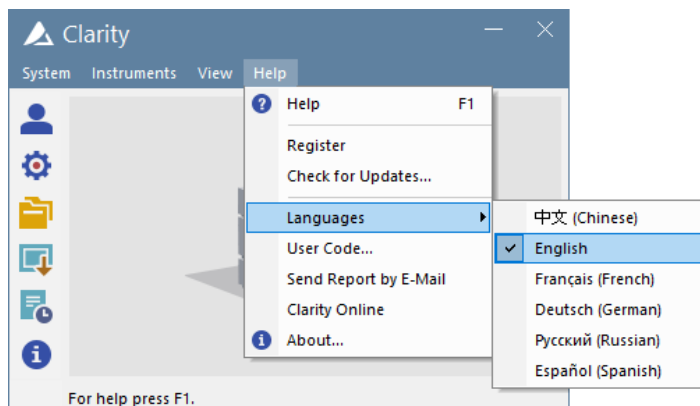


### Existing users

Registered Users of **Clarity (Lite) in English** can update Clarity to the latest version from our web page [www.dataapex.com](http://www.dataapex.com). This update is free of charge. Registration can be performed also at [www.dataapex.com](http://www.dataapex.com) or directly from Clarity.


There is no need to change neither a HW key (dongle) nor user code. Only if the original version is older than 2.4 a HW key has to be ordered (p/nC600).

After the update to the latest version a user can select the language directly from the SW:

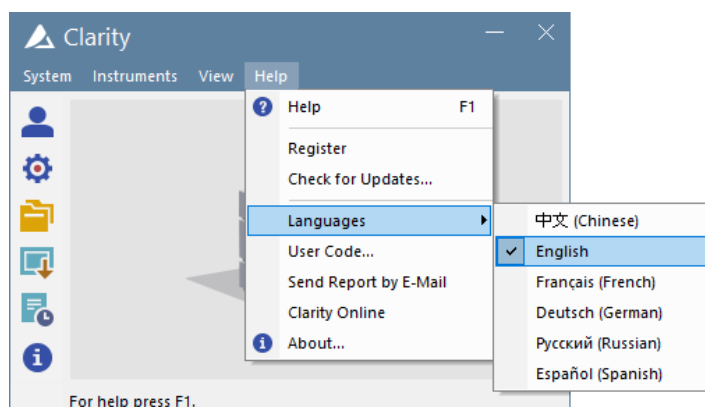

## How to upgrade from Chinese Clarity (p/n C50-CHS) or Clarity Lite (p/n C40-CHS) to Multilingual version

Users of **Clarity (Lite) in Chinese** need to purchase an upgrade (p/n C560 or C565) to Multilingual Clarity (Lite) Chromatography Software. **Clarity in Chinese** uses a different U/C which enables Chinese language only. A new standard user code issued for an upgrade will enable to use other languages

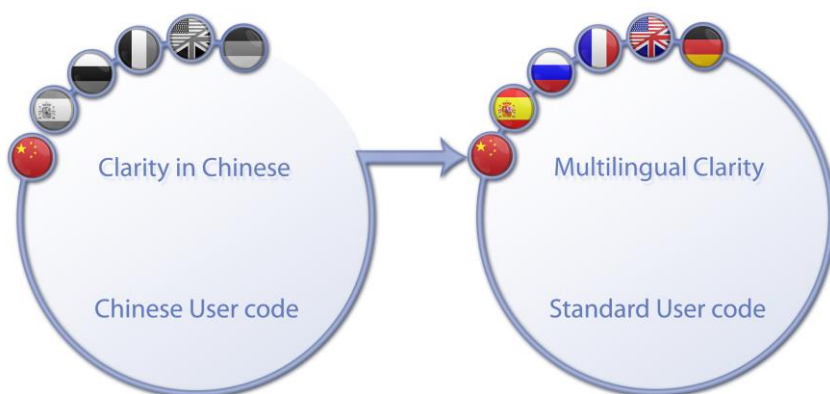

*There is also an upgrade path for users of obsolete chromatography SW CSW/Apex to Clarity (Lite) Chromatography SW. Users that will purchase an upgrade will automatically get the latest version of multilingual Clarity.*

## How to recognize user codes by the initial letter

<b>C</b>	Standard User Code	<b>Multilingual Clarity</b>
<b>L</b>	Clarity Lite User Code	<b>Multilingual Clarity Lite</b>
<b>X</b>	Chinese User Code	<b>Clarity in Chinese only</b>
<b>W</b>	Chinese Clarity Lite User Code	<b>Clarity Lite in Chinese only</b>