

## Clarity2Go

Analysis monitoring for smartphones

Clarity2Go is a mobile companion to Clarity Chromatography Station.

It enables to monitor selected parameters of running sequences or single analyses conveniently from your mobile. Its use is **free** for all customers with Clarity station.

*The monitoring works only if both (smartphone and PC with Clarity) have access to the internet.*

The whole solution consists of three parts:

- Clarity Software (Full, minimal version 4.0.4) communicating with the web server.
- Clarity2Go web server run by DataApex is provided to Clarity2Go users free of charge.
- Clarity2Go application for smartphones and tablets.

Clarity2Go application is currently available for Android ([Google Play](#)) and iOS ([Apple Store](#)).


For more information go to <http://www.dataapex.com/clarity2go>



## How to use Clarity2Go

### Enabling Clarity2Go in Clarity

In Clarity software on your PC use the *System - Clarity2Go...* command to invoke the *Clarity2Go Configuration* dialog.

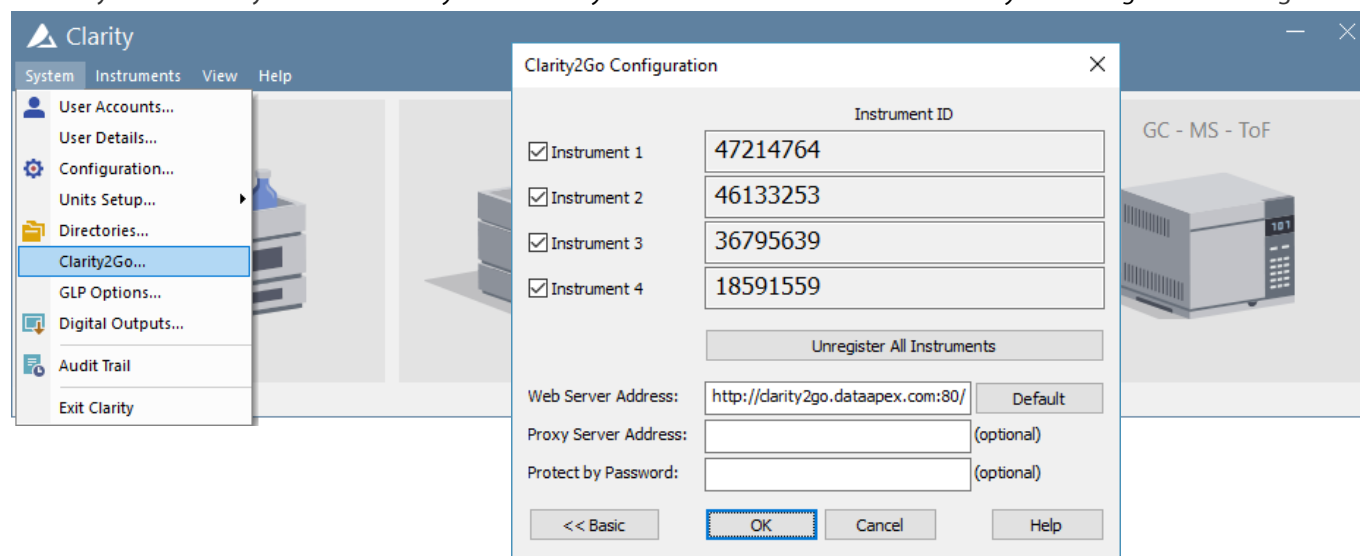

Clarity2Go Configuration	
Instrument ID	
<input checked="" type="checkbox"/> Instrument 1	47214764
<input checked="" type="checkbox"/> Instrument 2	46133253
<input checked="" type="checkbox"/> Instrument 3	36795639
<input checked="" type="checkbox"/> Instrument 4	18591559
Unregister All Instruments	
Web Server Address:	http://clarity2go.dataapex.com:80/ Default
Proxy Server Address:	(optional)
Protect by Password:	(optional)
<< Basic OK Cancel Help	

Select the instruments for monitoring, in the above example we have selected all 4 instruments.

After checking the checkbox the Instrument is automatically registered on our server and you will receive unique *Instrument ID* for it. Use these *Instrument IDs* in Clarity2Go application on your smartphone/tablet.

### Clarity2Go on your smartphone and tablet

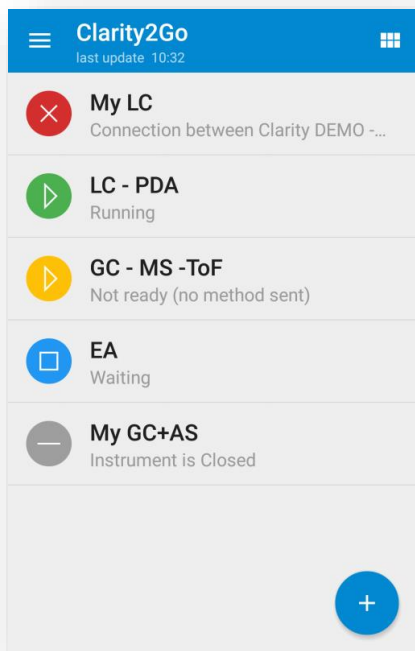


Figure 1-2: Android smartphone

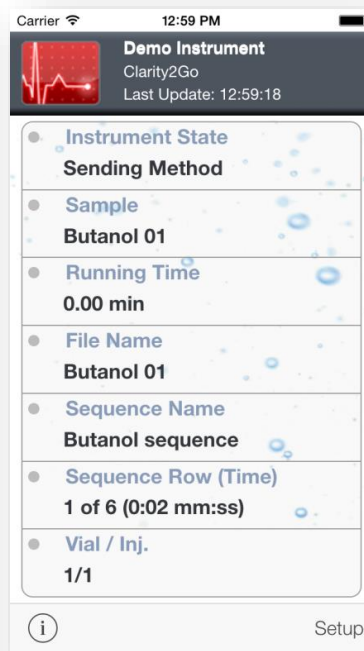
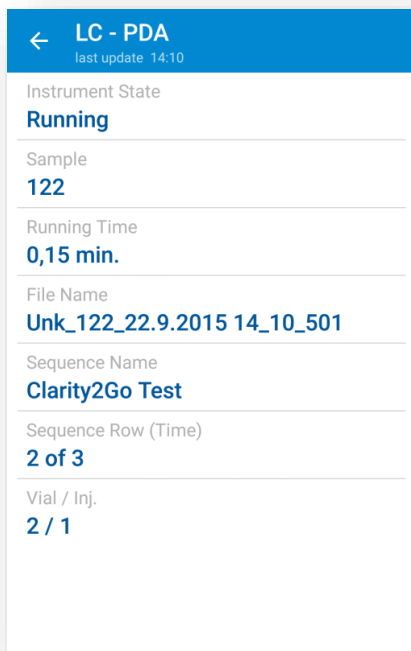


Figure 3: iOS smartphone

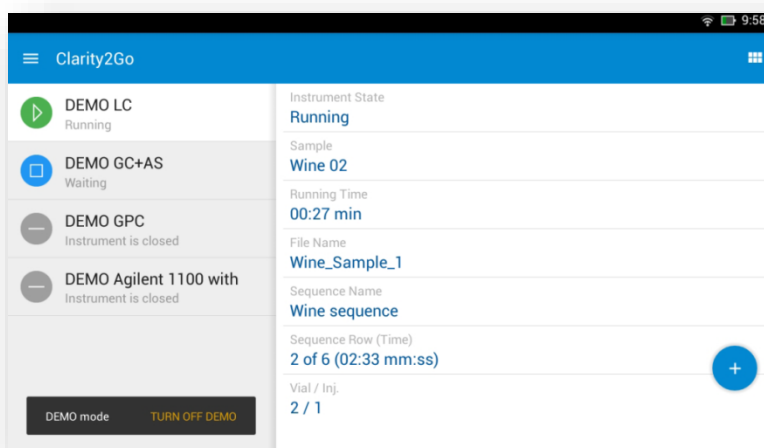


Figure 4: Android tablet

Clarity2Go was released with only few features and further development will follow the needs of Clarity users. If you have any suggestions or encounter any inconvenience, let us know at [support@dataapex.com](mailto:support@dataapex.com).

Help us make Clarity2Go better and give us your feedback!